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UviLine Series

Gamme Uviline

SPECTROPHOTOMETERS / SPECTROPHOTOMETRE

UviLine Series

The most current version of the present operating manual is available on our web site:

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<http://www.aqualabo-group.com/>

REMARKS:

The information contained in this document may be the object of patents or patent applications by Aqualabo Analyse. The possession of this document does not confer any license to such patents.

The following names are Aqualabo Analyse's registered trademarks throughout the world:

UviLine

This manual is updated periodically. The updates are included in the new editions.

All information supplied in this edition of the manual may be amended before the products described herein are available.

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Version française : page 117 to page 207

Copyright ©

Ales 2016, Aqualabo Analyse
Reproduction in whole - or even in part - is prohibited without the express written permission of Aqualabo.

CONTENTS

1	WARRANTY	9
2	INFORMATION.....	11
3	SAFETY.....	13
3.1	SAFETY INFORMATION	13
3.1.1	SAFETY INFORMATION IN THE OPERATING MANUAL.....	13
3.1.2	SAFETY SIGNS ON THE PRODUCT	13
3.2	SAFE OPERATION	13
3.2.1	AUTHORIZED USE.....	13
3.2.2	REQUIREMENTS FOR SAFE OPERATION.....	13
3.2.3	UNAUTHORIZED USE	13
3.3	HANDLING OF HAZARDOUS SUBSTANCES.....	14
3.4	GENERAL NOTES ON HANDLING.....	14
3.5	PACKING.....	14
4	OVERVIEW.....	15
4.1	FRONT OF THE INSTRUMENT	15
4.2	SOCKET FIELD ON THE REAR PANEL	15
4.2.1	USB-A PORT	16
4.2.2	USB-B PORT	16
4.3	KEYBOARD.....	17
4.3.1	OVERVIEW	17
4.3.2	KEY FUNCTIONS.....	17
4.3.3	FUNCTION KEYS.....	18
4.3.4	ALPHANUMERIC KEYBOARD	18
4.3.4.1	Characters sets	18
4.3.4.2	Command principle.....	18
4.3.4.3	Special characters	18
4.3.4.4	How to correct a bad character.....	18
4.4	DISPLAY.....	19
4.4.1	DISPLAY ELEMENTS	19
4.4.2	SYMBOLS IN THE STATUS LINE	19
4.5	CUVETTE COMPARTMENT.....	20
5	TECHNICAL DATA	21
5.1	MEASUREMENT CHARACTERISTICS.....	21
5.2	MEASUREMENT MODE	24
5.2.1	CONCENTRATION.....	24
5.2.2	ABSORBANCE / % TRANSMISSION.....	24
5.2.3	MULTI WAVELENGTHS	24
5.2.4	SPECTRUM.....	24
5.2.5	KINETICS.....	24
5.3	MEMORY FOR MEASURED VALUES	25
5.3.1	MEMORY CAPACITY.....	25
5.3.2	OUTPUT OPTIONS	25
5.3.3	FILE FORMATS	25
5.4	GENERAL INSTRUMENT DATA	25
5.5	AVAILABLES LANGUAGES:	26
6	COMMISSIONING	27
6.1	BUFFER BATTERIES INSTALLATION	27
6.2	CONNEXION OF POWER SUPPLY	27
6.3	CHOICE OF THE LANGUAGE	28
6.4	DATE AND TIME SETTING UP	28

UviLine Series

6.4.1	SELF-TEST.....	29
6.4.2	WARM-UP TIME.....	29
7	OPERATION.....	31
7.1	SWITCHING ON OR OFF THE PHOTOMETER	31
7.2	GENERAL OPERATING PRINCIPLES.....	32
7.2.1	NAVIGATING	32
7.2.2	CUSTOMER METHODS AND PROFILE IMPORTATION/EXPORTATION.....	33
7.2.3	PDF FILE PRINTING.....	33
8	CONCENTRATION MODE	35
8.1	GENERAL POINT.....	35
8.2	METHOD PROGRAMMING	35
8.3	METHOD SAVING.....	38
8.4	METHOD DELETING	39
8.5	METHOD EXECUTION	39
8.5.1	WITH STANDARD SOLUTIONS.....	39
8.5.1.1	Measure.....	39
8.5.1.2	Calibration curve displaying.....	41
8.5.1.3	Point deleting.....	42
8.5.1.4	Calibration curve printing.....	43
8.5.1.5	Sample measurement	43
8.5.1.6	Results printing.....	44
8.5.1.7	Sample dilution.....	44
8.5.1.8	Sample blank value	45
8.5.2	WITH VALUE PAIRS.....	46
8.5.3	WITH A FORMULA.....	47
8.6	METHOD EDITING.....	47
9	ABSORPTION / TRANSMISSION MODE.....	49
9.1	GENERAL POINTS	49
9.2	MEASUREMENT.....	49
9.3	RESULTS PRINTING	49
10	SPECIAL/ MULTI-WAVELENGTHS.....	51
10.1	BASIC INFORMATION.....	51
10.2	METHOD PROGRAMMING	51
10.3	METHOD DELETING	53
10.4	METHOD SELECTION.....	53
10.5	METHOD EXECUTION	54
11	SPECTRUM MODE.....	56
11.1	GENERAL POINTS	56
11.2	METHOD PROGRAMMING	56
11.3	METHOD SAVING.....	57
11.4	METHOD DELETING	57
11.5	METHOD SELECTION.....	57
11.6	MEASUREMENT EXECUTION.....	57
11.7	SPECTRA TREATMENT	58
11.7.1	DOWNLOADING OF STORED SPECTRUM.....	58
11.7.2	CURSOR.....	58
11.7.3	ZOOM.....	58
11.7.4	MANUAL CHANGE OF ABSORBANCE SCALE.....	60
11.7.5	CALCULATIONS.....	60
11.7.5.1	Picks and valleys detection.....	60
11.7.5.2	Mark points.....	60
11.7.5.3	Deleting of mark points.....	61
11.7.5.4	Area calculation	62
11.7.5.5	Derivative calculation.....	62
11.7.5.6	Spectra comparison.....	63

11.7.5.7	Addition of two spectra	63
11.7.5.8	Subtraction of one spectrum from another one.....	64
11.7.5.9	Division of a spectrum	64
11.7.5.10	Addition of a constant	64
11.7.5.11	Multiplication by a constant.....	65
12	KINETICS MODE.....	67
12.1	GENERAL POINTS	67
12.2	METHOD PROGRAMMING	67
12.3	METHOD SELECTION	68
12.4	METHOD SAVING	68
12.5	METHOD DELETING	69
12.6	LIMITATION OF METHODS LIST	69
12.7	METHOD EXECUTION	70
12.8	METHOD EDITING	71
12.9	KINETICS TREATMENT	71
12.9.1	<i>KINETIC CURVE PRINTING</i>	71
12.9.2	<i>CURSOR</i>	71
12.9.3	<i>KINETIC CURVE SAVING</i>	71
12.9.4	<i>DOWNLOADING OF STORED KINETIC CURVE</i>	71
12.9.5	<i>« EDIT » MENU</i>	72
12.9.5.1	Slope	72
12.9.5.2	Comparison of kinetic curves.....	72
12.9.5.3	Subtraction of a kinetic curve from another one	73
12.9.6	<i>MANUAL CHANGE OF ABSORBANCE SCALE</i>	73
12.10	CURVE PRINTING	73
13	MEMORY.....	75
13.1	OVERVIEW	75
13.2	INSTRUCTIONS ON USING USB MEMORY DEVICES	76
13.3	MEASUREMENT DATASETS	76
13.3.1	<i>ELEMENT OF A MEASUREMENT DATASET</i>	76
13.3.2	<i>OPERATIONS WITH MEASUREMENT DATASETS</i>	76
13.3.3	<i>RESULT RECORDING</i>	77
13.3.3.1	Manual recording.....	77
13.3.3.2	Automatic recording with « AutoStore » function	77
13.3.4	<i>RESULT DISPLAYING</i>	78
13.3.5	<i>RESULT FILTERING</i>	80
13.3.5.1	Filtering of a results group	80
13.3.5.2	Filters inversion	80
13.3.5.3	Erase stored measurement dataset.....	80
13.3.5.4	Result deleting.....	81
13.3.5.5	Saving Kinetic recordings, spectrUM.....	81
13.3.5.6	Saving data as a pdf file	81
13.3.6	<i>SAVING / EXPORTING FILES</i>	81
13.3.6.1	Copying all measurement data files to a USB memory device	81
13.3.6.2	Copying user-defined methods / profiles to a USB memory device.....	82
13.3.6.3	Copy of individual file on USB memory	82
13.3.7	<i>PRINTING THE DATA (USB)</i>	83
13.3.8	<i>SETTING FOR DATA TRANSMISSION</i>	83
13.3.9	<i>IMPORTING FILES</i>	84
13.3.9.1	Importing spectra or kinetic recordings from a USB memory device	84
13.3.9.2	Importing methods / profiles from a USB memory device.....	84
14	USERS MANAGEMENT	85
14.1	USER LEVEL AND USER RIGHTS	85
14.2	ACTIVATION/DEACTIVATION OF USER MANAGEMENT	85
14.2.1	<i>ACTIVATION OF USER MANAGEMENT FUNCTION</i>	86
14.2.2	<i>DEACTIVATION OF USER MANAGEMENT FUNCTION</i>	86
14.3	CREATION, MODIFICATION OR DELETING A USER ACCOUNT	86
14.3.1	<i>USER ACCOUNT CREATION</i>	86

UviLine Series

14.3.2	USER ACCOUNT MODIFICATION	87
14.3.3	PASSWORD MODIFICATION	87
14.3.4	USER ACCOUNT SUPPRESSION	87
15	VARIOUS ADJUSTMENT	105
15.1	LANGUAGE CHOICE	105
15.2	DATE AND TIME ADJUSTMENT	105
15.3	CONTRAST ADJUSTMENT	105
15.4	RESET	106
15.5	LAMP SERVICING COUNT DOWN	106
15.6	SPECTROPHOTOMETER INFORMATION ([INFO])	106
15.7	TIMER	107
15.8	SOFTWARE UPDATING	107
16	MAINTENANCE.....	109
16.1	LAMP CHANGING	109
16.2	BUFFER BATTERIES CHANGING	110
16.3	CLEANING	111
16.3.1	CLEANING YOUR SEPCTROPHOTOMETER.....	111
16.3.2	CLEANING THE LENSES	111
16.4	WHAT CAN YOU DO, IF.....	111
16.4.1	WARNING IF CUVETTE IS BROKEN.....	111
16.5	TROUBLESHOOTING	112
17	ACCESSORIES	113
17.1	10MM CELL HOLDER (REF. 70VI0604).....	113
17.2	PORTE CUVE JUSQU'A 100MM (REF. 70VI0605)	113
17.3	PORTE CUVE 10MM THERMOSTATABLE (REF. 70VI0602)	113
17.4	PORTE TUBE 16MM (REF. 70VI0607)	113
17.5	AUTOMATIC CELL CHANGER TURRET (REF. 70VI0600).....	113
17.6	SIPPER (REF. 70VI0601).....	114
17.7	EFFET PELTIER (REF. 70VI0603)	114
17.8	SIPPER WITH PELTIER EFFECT (REF. 70VI0610)	114
17.9	DIAPHRAGMS FOR MICRO-CELLS (REF. 70VI0608)	115

1 WARRANTY

UviLine Series instruments sold by Aqualabo Analyse is guaranteed against any manufacturing defects for three years (unless otherwise stated by AQUALABO ANALYSE) with effect:

- From the technical acceptance of the equipment in the factory by the buyer or his designee,
- or failing this :
 - * For Metropolitan France: from the date on the delivery note.
 - * For other destinations: from the date of factory shipment certified by air waybill, consignment note or bill of lading.

Our company guarantee applies exclusively to defectiveness arising from a design fault or from a concealed defect. It is strictly limited to the free dispatching of replacement parts (except for consumable items) or to the repairing of the equipment in our workshops within a deadline of 10 working days (shipping delay not included).

By express agreement, the following are strictly excluded from our guarantee:

- All damages, notably for staff costs, loss of earnings, business trouble, etc
- Any breakdown due to an incorrect use of the equipment (non adapted mains, fall, attempt at transformation, etc) or to a lack of maintenance by the user or to poor storage conditions.
- Any breakdown due to the use of parts not supplied by AQUALABO ANALYSE, on AQUALABO ANALYSE equipment
- Any breakdown due to the transporting of the equipment in packaging which is not its original packaging
- The lamps, the cells and generally any item which appears in the "accessories" section on the price list.

Our customers are kindly asked to apply for our consent before returning any instrument for repair. No return of materials may be accepted without the prior written consent of our Servicing after Sales Management which will precise the terms of such return. If the above consent is given, articles shall be returned in their original packaging on a prepaid basis to the following address:

AQUALABO ANALYSE – USINE SECOMAM - 91 avenue des Pins d'Alep 30100 ALES - France

We reserve the right to reship all instruments received collect failing such consent.

Whatever method and conditions of transport are chosen for the shipment of the equipment to be repaired under guarantee, in the original packaging, the corresponding costs and the insurance costs will be payable by the customer.

Any damage connected to the return transport of the equipment falls within the framework of the guarantee on the express condition that the customer has sent his complaint within forty-eight hours by registered letter with acknowledgement of receipt to the carrier. A copy of the letter should be sent to AQUALABO ANALYSE.

For equipment with a guaranty card, this is only applicable if the card delivered with the equipment is returned to AQUALABO ANALYSE duly completed.

SOFTWARE GUARANTEE

The software is guaranteed by the designer or the distributor of the software under the conditions specified in the literature accompanying the aforementioned software packages.

Under no circumstances whatsoever will AQUALABO ANALYSE supply any type of guarantee for software packages.

By express agreement, all damages, notably for staff costs, lost of earnings; business trouble, etc are strictly excluded from our guarantee.

The customer is informed that the software cannot be guaranteed exempt from defects or bugs.

TRADE SECRET AND PROPERTY RIGHTS

This document is protected by AQUALABO ANALYSE copyright (c) 2003 and the copying rights are explicitly reserved. The software supplied with the equipment or referenced contains trade secrets and confidential information which are AQUALABO ANALYSE's property. It is legally protected by the international copyright (c) laws.

AQUALABO ANALYSE grants a license to use its software to the user. This may not be disclosed, used or duplicated with the intention to save it, without AQUALABO ANALYSE's written permission. The beneficiary must attach a copy of this document to all authorized partial or total reproductions.

2 INFORMATION

The AQUALABO ANALYSE equipment has been designed, manufactured, tested and inspected according to the ISO 9001 standards (Version 2000).

If the unit is not immediately installed, it should be stored in a dry and clean area. The storage temperature should be between 10 and 35°C.

AQUALABO ANALYSE equipment is carefully inspected before it is packed. As soon as you receive your equipment, check the condition of the packaging and if you notice any problems, notify your carrier within 48 hours. Then consult the packing list and check that everything is in order. Finally, if you discover that something is missing, or if the goods are damaged immediately notify AQUALABO ANALYSE.

The main differences between UviLine 9300 and UviLine 9600 spectrophotometers are the wavelengths ranges and the light sources.

UviLine 8100C / 9100C / 9300

Wavelengths range: 320 – 1100 nm
Light source: Halogen lamp 5VDC 10W

UviLine 9400C / 9600

Wavelengths range: 190 – 1100 nm
Light source: Xenon lamp.

UviLine Series instruments are fully designed and manufactured by AQUALABO ANALYSE in Ales France.

AQUALABO ANALYSE / SECOMAM

91 Avenue Des Pins D'Alep • 30100 Alès • France

Tel : +33 (0)4 66 54.26.95

Fax : +33 (0)9 72 35 16 80

E-mail : info@aqualabo.fr

Web: www.aqualabo-group.com

3 SAFETY

1.1 SAFETY INFORMATION

1.1.1 SAFETY INFORMATION IN THE OPERATING MANUAL

This operating manual provides important information on the safe operation of the product. Read this operating manual thoroughly and make yourself familiar with the product before putting it into operation or working with it. The operating manual must be kept in the vicinity of the product so you can always find the information you need.

Important safety instructions are highlighted in this operating manual. They are indicated by the warning symbol (triangle) in the left column. The signal word (e.g. "CAUTION") indicates the level of danger:

WARNING

Indicates a possibly dangerous situation that can lead to serious (irreversible) injury or death if the safety instruction is not followed.

1.1.2 SAFETY SIGNS ON THE PRODUCT

Note all labels, information signs and safety symbols on the product. A warning symbol (triangle) without text refers to safety information in this operating manual.

1.2 SAFE OPERATION

1.2.1 AUTHORIZED USE

The authorized use of the photometer consists exclusively of the carrying out of photometric measurements according to this operating manual. Follow the technical specifications of the cells. Any other use is considered **unauthorized**.

1.2.2 REQUIREMENTS FOR SAFE OPERATION

Note the following points for safe operation:

- The product may only be operated according to the authorized use specified above.
- The product may only be supplied with power by the energy sources mentioned in this operating manual.
- The product may only be operated under the environmental conditions mentioned in this operating manual.
- The product may not be opened.

1.2.3 UNAUTHORIZED USE

The product must not be put into operation if:

- it is visibly damaged (e.g. after being transported)
- it was stored under adverse conditions for a lengthy period of time

User qualification

Carrying out photometric determinations with the aid of test sets frequently requires the handling of hazardous substances.

We assume that the operating personnel know how to handle hazardous substances due to their professional training and experience. The operating personnel must particularly be able to understand and correctly implement the safety labels and safety instructions on the packages and inserts of the test sets.

1.3 HANDLING OF HAZARDOUS SUBSTANCES

When developing test sets, Orchidis carefully sees that the tests can be carried out as safely as possible. Some hazards by dangerous substances, however, cannot always be avoided.

If self-produced tests or solutions are used, the responsibility concerning any risks caused by those tests or solutions lies with the user (personal responsibility).

WARNING

Improper handling of certain reagents can cause damage to your health. In any case follow the safety labels on the packing and the safety instructions of the package insert. Protective measures specified there have to be followed exactly.

Safety datasheets

The safety datasheets of the chemicals comprise all instructions on safe handling occurring hazards, preventive actions and actions to take in hazardous situations. Follow these instructions in order to work safely.

1.4 GENERAL NOTES ON HANDLING

The UviLine instruments are an optical precision meters. Therefore, it should always be handled with care, especially in mobile use.

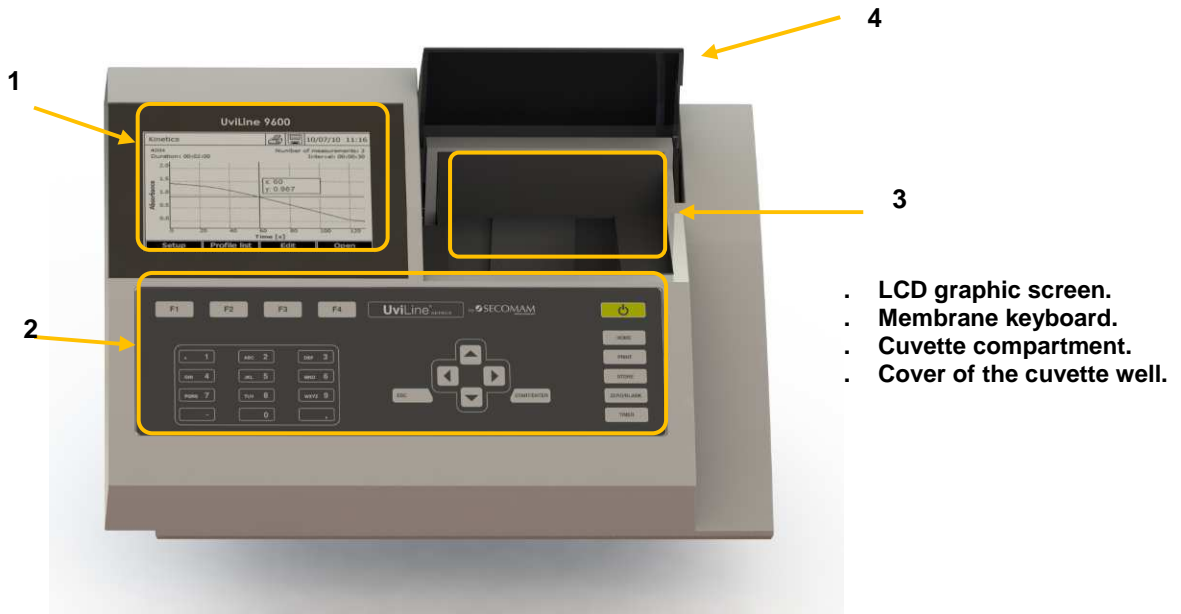
- Always protect the meter from conditions that could damage the mechanical, optical and electronic components. Heed the following points especially:
- The temperature and humidity during operation and storage must be within the limits specified in chapter TECHNICAL DATA.
- The following influences always have to be avoided with the meter:
 - Extreme dust, moisture and wetness
 - Exposure to intensive light and heat
 - Fumes that are corrosive or contain high concentrations of solvents.
- For measuring, the meter must be placed on a flat surface.
- Spilled liquid or other material should be removed immediately
- If a cell has broken in the cell shaft, the cell shaft should be cleaned immediately
- The cell shaft should always be closed when the photometer is not used.
- For the instrument to be transported the cell shaft has to be empty.
- For mobile used we recommend using the suitable transport case
- Always make sure that the instrument is connected on the good voltage. (Between 100 – 240V 50-60Hz)
- When dangerous substances for health and environment are used, the laboratory or site rules, where the instrument is installed must be followed.
- Take all the necessary precautions, during the use the instrument, to protect the operator from eventual liquids leaks or spills or possible radiations (protective gloves, anti-UV radiation glasses , protected clothes, etc)
- Keep the sample compartment clean.
- The Xenon lamp used in the UviLine 9400C & 9600 emits UV radiation.
- Install the instrument in a ventilated area because it is likely to generate ozone, which, beyond the limits below, can harm health.
 - Exposure average value = 100 ppb
 - Exposure limits value = 200 ppb
- All operations made inside the instrument, must be done by AQUALABO ANALYSE or by AQUALABO ANALYSE's authorized technicians.

1.5 PACKING


The spectrophotometer is sent out in a protective transport packing. Keep the original packing including the inner packing to protect the instrument against hard shocks if it has to be transported. The original packing is also required for the proper return of the instrument if it has to be repaired. Note that damage caused by improper transport voids all warranty claims.

4 OVERVIEW

1.6 FRONT OF THE INSTRUMENT

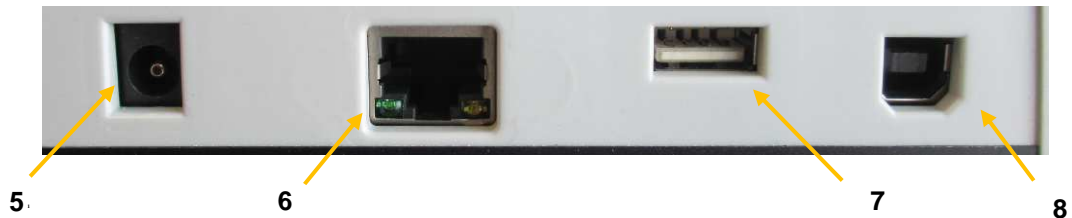1.7 SOCKET FIELD ON THE REAR PANEL

You can connect the following accessories to the photometer:

- PC (USB-B)
- Printer (USB-A → The data is printed out with <PRINT>)
- USB memory device (USB-A)
- USB-PC keyboard (USB-A)

- 5. Connection for power supply.
- 6. Ethernet RJ45.
- 7. USB-A connection.
- 8. USB-B connection.

Suitable are all printers that can interpret the PCL-3 printer control language.

If you want to connect several USB devices such as a USB-PC keyboard and a USB memory device to the meter, you can increase the number of USB-A sockets by a commercially available USB-2 hub with separate power supply.

1.7.1 USB-A PORT

The USB-A port allows:

- To print data when a printer is connected. With <PRINT> key data are sent on the printer.
- To records data on USB memory (USB key) when it is connected.
- To update spectrophotometer software and methods data.
- To connect an external USB keyboard.

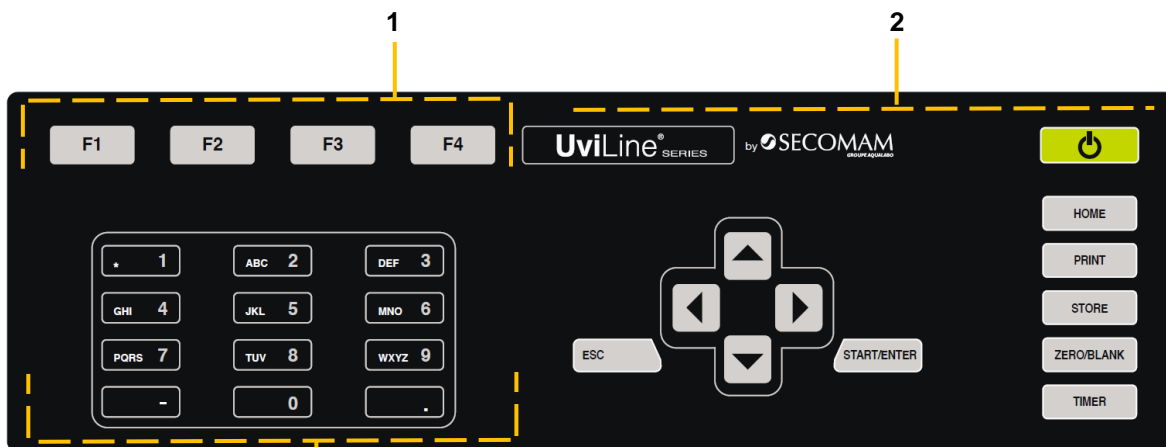
1.7.2 USB-B PORT

Allows the spectrophotometer and PC to be connected directly by emulating a RS232 port. A USB ↔ External RS232 adapter is not required to connect the spectrophotometer to a free PC RS232C port. Please contact Orchidid Aqualabo Analyse for the driver package.

UviLine Series

1.8 KEYBOARD

1.8.1 OVERVIEW

1. Function keys F1 to F4 (functions menu depending)
2. Fixed key functions.
3. Alphanumeric keys block.

1.8.2 KEY FUNCTIONS

Key	Designation	Function
	<ON/OFF>	Switch the spectrophotometer ON and OFF
	<HOME>	Switches to the main menu from any operating situation. Actions that are not completed are cancelled.
	<PRINT>	Downloads the displayed value to an interface.
	<STORE>	Saves a displayed value or spectrum or kinetic curve.
	<ZERO BLANK>	Starts one of the following measurements, depending on the operating situation: - Zero adjustment - Blank value measurement - Baseline measurement
	<TIMER>	Open "Timer" menu.
	<ESC>	Cancels the running action. - Entries that have not yet been accepted are discarded. - Switches to the next higher menu level.
	<START ENTER	- Starts an action (e.g. measurement) - Opens a selected menu - Confirms a selection or entry
	<▲> or <▼>. <◀>	Moves the selection in menus and lists one position up or down. - Deletes the character left of the cursor during character entries - Moves the cursor to the left in a spectrum or kinetic diagram.

– Moves the cursor to the right in a spectrum or kinetic diagram.

1.8.3 FUNCTION KEYS

The function keys F1 to F4 have different functions depending on the operating situation. The current functions are displayed in the function key menu at the bottom edge of the display.

1.8.4 ALPHANUMERIC KEYBOARD

Numerals, letters, punctuation marks and special characters are entered with the alphanumeric keypad of the meter or using an external keyboard.

1.8.4.1 Characters sets

The following characters are available:

- Numerals (0 ... 9)
- Letters (A ... Z) and (a ... z).
- Punctuation marks (. and -)
- Special characters ° / + Δ, Σ, μ, #, %, : et ()

1.8.4.2 Command principle

Entering characters is always possible if there is an input field on the display.

The numerals and characters (except for the small letters) assigned to the keys of the alphanumeric keypad are printed on the keys.

Example: With the <2/ABC> key you can enter the following characters: 2, A, B, C, a, b, c.

Select the required character by pressing the key several times (similar to a mobile phone). When pressing a key that is assigned to several characters once, the respective numeral appears first.

To enter a numeral, one key pressing is always sufficient.

When pressing the key for the first time a line pops up that displays all characters possible with this key. The currently selected character is highlighted.

A character is taken over in the input field if

- the character is highlighted for more than one second,
- the character is confirmed with <START-ENTER>,
- Another alphanumeric key is pressed.

Remark

During mere number entries (such as entering a wavelength), the keys of the alphanumeric keypad are assigned to the respective numeral only. Each pressing key directly enters the numeral (like a pocket calculator).

1.8.4.3 Special characters

To enter special characters, use <1/*> key.

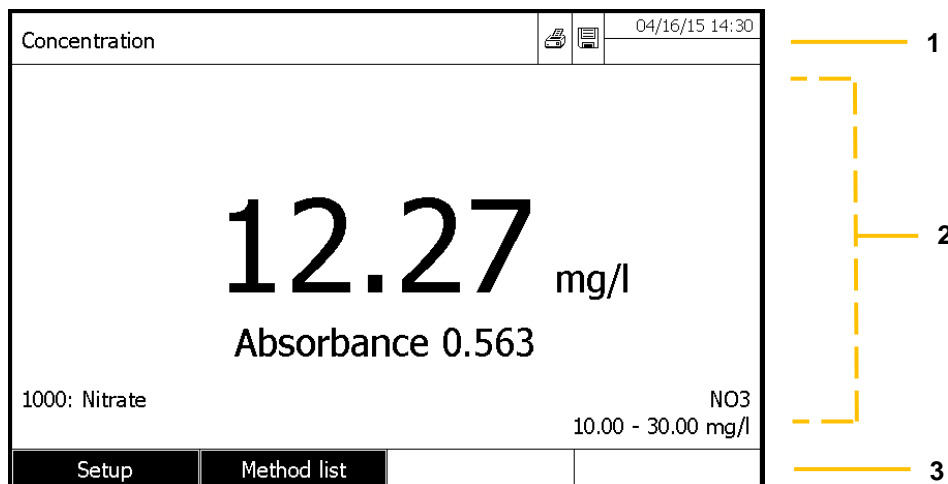
1.8.4.4 How to correct a bad character

Use <◀> key, to delete all characters until THE INCORRECT one and start again the entries from there.

UviLine Series

1.9 DISPLAY

1.9.1 DISPLAY ELEMENTS


1. Status line (current state, date and time)
2. Display range for menus and measurement results
3. Function keys menu

1.9.2 SYMBOLS IN THE STATUS LINE

Symbol	Designation	Function
	Save	The <STORE> key is active. You can store the displayed data with <STORE>
	Printer	The <PRINT> key is active. You can output to an interface the displayed data with <PRINT>


1.10 CUVETTE COMPARTMENT

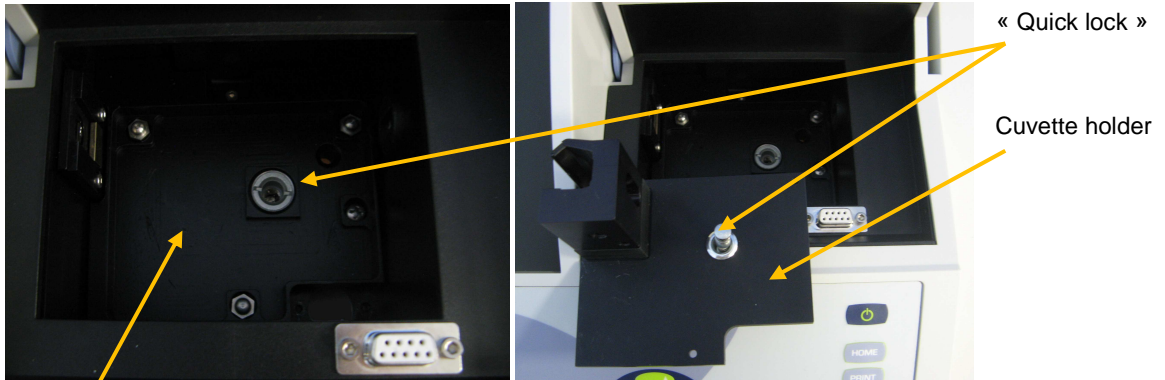The cuvette holder is quickly install thank to a special system named "Quick lock"

Cuvette compartment

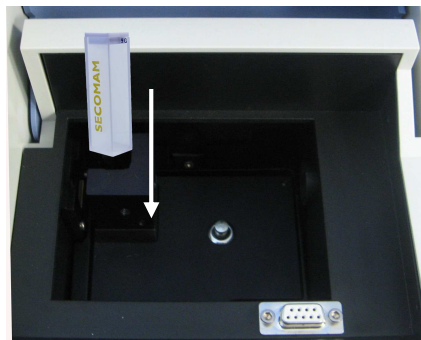- Set in position the cuvette holder (Quick lock system above the hole inside the cuvette well).
- Push, with finger, the "quick lock" system.
- Release the "quick lock" system.
- The cuvette holder is installed. And ready to receive a cuvette
- The beam is going from the left to the right.

5 TECHNICAL DATA

1.11 MEASUREMENT CHARACTERISTICS

Uviline 8100C

Wavelength range	• 320-1100 nm
Light Source	• Halogen
Bandpass	• 6 nm
Incremental WL step	• Reading: 0.1 nm • Setting : 1 nm
Wavelength accuracy	• ± 2 nm
Wavelength repeatability	• $\pm 0,5$ nm
Absorbance range	• $\pm 3,300$
Absorbance resolution	• 0,001
Photometric accuracy	• $\pm 0,005$ Abs (0,5 Abs) • $\pm 0,008$ Abs (1,0 Abs) • $\pm 0,016$ Abs (2,0 Abs)
Stray light	• $<0.1\%$ at 340nm (NaNO ₂)
Flatness baseline	• $\pm 0,010$ Abs
Scanning speed	• Low – medium – fast
Update	• Via USB port
Interface	• 1 USB-A, 1 USB-B, 1 RS 232C
IP standard	• IP 30 with drain in the cell compartment
Power supply	• 110-220V 50/60 Hz - specific country cable(s)
Dimensions (L x W x H)	• 404 x 314 x 197 mm
Weight (net)	• 4,7 Kg
Warranty	• 3 years


	UviLine 9100 Visible unit	UviLine 9400 UV-Visible unit
Wavelength range	• 320-1100 nm	• 190-1100 nm
Light Source	• Halogen	• Xenon
Bandpass	• 4 nm	• 4 nm
Incremental WL step	• Reading: 0.1 nm • Setting : 1nm	• Reading: 0.1 nm • Setting : 1nm
Wavelength accuracy	• ± 1 nm	• ± 1 nm
Wavelength repeatability	• ± 0,2 nm	• ± 0,2 nm
Absorbance range	• ± 3,300	• ± 3,300
Absorbance resolution	• 0,001 Abs ou 0,1% T	• 0,001 Abs ou 0,1% T
Photometric accuracy	• ± 0,003 Abs (0,5 Abs) • ± 0,005 Abs (1,0 Abs) • ± 0,010 Abs (2,0 Abs)	• ± 0,003 Abs (0,5 Abs) • ± 0,005 Abs (1,0 Abs) • ± 0,010 Abs (2,0 Abs)
Stray light	• ≤ 0.1 % to 340 nm (GG375) 400 nm (GG408)	• ≤ 0.1 % to 220 nm (NaI) 340 nm (GG375) 400 nm (GG408)
Flatness baseline	• ± 0,010 Abs	• ± 0,010 Abs
Scanning speed	• Low – medium – fast	• Low – medium – fast
Update	• Via USB port	• Via USB port
Interface	• 1 USB-A, 1 USB-B, 1 RS 232C	• 1 USB-A, 1 USB-B, 1 RS 232C
IP standard	• IP 30 with drain in the cell compartment	• IP 30 with drain in the cell compartment
Power supply	• 110-220 V 50/60 Hz specific country cable	• 110-220 V 50/60 Hz specific country cable
Température (°C) / Humidity	• Use: 10 to 35 °C • Storage -25 °C to 65 °C	• Use: 10 to 35 °C • Storage -25 °C to 65 °C
Dimensions (L x W x H)	• 404 x 197 x 314 mm	• 404 x 197 x 314 mm
Weight (net)	• 4 Kg	• 4 Kg
Warranty	• 2 years	• 2 years

UviLine Series

	9300	9600
Wavelengths range	• 320-1100 nm	• 190-1100 nm
Light source	• Halogen	• Xenon
Spectral bandwidth	• 4 nm	• 4 nm
Incremental WL step	• Reading: 0.1 nm • Setting: 1 nm	• Reading: 0.1 nm • Setting: 1 nm
Wavelength accuracy	• ± 1 nm	• ± 1 nm
Wavelength repeatability	• $\pm 0,5$ nm	• $\pm 0,5$ nm
Absorbance range	• $\pm 3,500$ Abs	• $\pm 3,500$ Abs
Absorbance resolution	• 0,001 Abs ou 0,1% T	• 0,001 Abs ou 0,1% T
Photometric accuracy	• $\pm 0,003$ Abs (0,5 Abs) • $\pm 0,005$ Abs (1,0 Abs) • $\pm 0,010$ Abs (2,0 Abs)	• $\pm 0,003$ Abs (0,5 Abs) • $\pm 0,005$ Abs (1,0 Abs) • $\pm 0,010$ Abs (2,0 Abs)
Stray light	• < 0,1 %T at 340 nm (GG375) • < 0,1 %T at 408 nm (GG408)	• < 1 %T at 198nm (KCl) • < 0,1 %T at 220nm (NaI) • < 0,05 %T at 340 nm (GG375) • < 0,05 %T at 408 nm (GG408)
Flatness baseline	• $\pm 0,002$ Abs	• $\pm 0,002$ Abs
Scanning speed	• Low – medium – fast > 600 nm/min	• Low – medium – fast > 800 nm/min
Update	• Via USB port	• Via USB port
Interface	• 1 USB-A, 1 USB-B, Ethernet RJ45	• 1 USB-A, 1 USB-B, Ethernet RJ45
IP	• IP 30 with drain in the cell compartment	• IP 30 with drain in the cell compartment
Power Supply	• 110-220 V 50/60 Hz specific country cable	• 110-220 V 50/60 Hz specific country cable
Temperature (°C)	• Use: 10°C to 35°C • Storage: 25°C to 65°C	• Use: 10°C to 35°C • Storage: 25°C to 65°C
Dimensions (L x l x h)	• 404 x 197 x 314 mm	• 404 x 197 x 314 mm
Weight (net)	• 4 Kg	• 4 Kg
Warranty	• 3 years	• 3 years

The specifications above are valid with the cuvette cover closed. To ensure accuracy and performances make measurements in low and stable light envirement (as dark as possible).

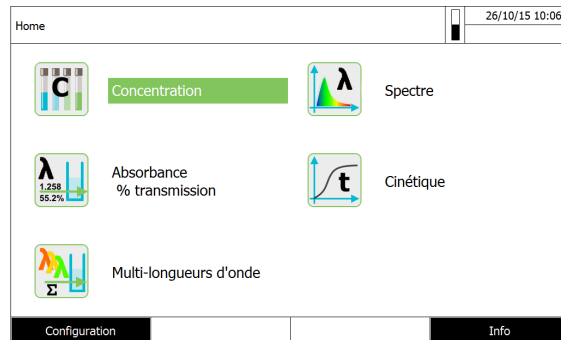1.12 MEASUREMENT MODE

1.12.1 CONCENTRATION

- Measurement with permanently programmed methods, adapted to the aqualabo test set program
- Program support for the creation of additional user-defined methods (max. 100)
- Citation forms and units method dependent
- Display of the absorbance value can be added

1.12.2 ABSORBANCE / % TRANSMISSION

- Measurement against own reference absorbance value possible

1.12.3 MULTI WAVELENGTHS

- Freely definable calculations from absorbance values at up to 10 different wavelengths
- Calculations can be stored as methods (max. 499)

1.12.4 SPECTRUM

- Absorbance or % transmission mode
- Limits freely selectable within the wavelength range
- Increment: 1 nm
- Settings can be stored as profiles (max. 20)
- Evaluation functions: Cursor scanning, zoom, min./max.recognition, peak area determination, derivation, smoothing, multiplication by constants, addition of constants, addition and subtraction of spectra, formation of the quotient of two spectra

1.12.5 KINETICS

- Absorbance or % transmission mode
- Minimal adjustable scan interval: 1 s (if the absorbance of the test sample is high, the scan interval is extended due to the longer duration of the individual measurements)
- Settings can be stored as profiles (max. 20)
- Evaluation functions: Cursor scanning, zoom, min/max determination, slope calculation (for an interval or total), enzymatic activity

1.13 MEMORY FOR MEASURED VALUES

1.13.1 MEMORY CAPACITY

- 5000 single measured values from the measuring modes, concentration, absorbance/ % transmission and multi wavelengths
- 40 MByte internal memory, sufficient for approx. 500 spectra and 400 kinetic curves (sample values based on the following assumptions: All spectra over a wavelength range of 600 nm and all kinetic curves with 150 single values each)

1.13.2 OUTPUT OPTIONS

USB memory device, printer, PC

1.13.3 FILE FORMATS

ASCII, *.csv

1.14 GENERAL INSTRUMENT DATA

Dimensions	404 x 197 x 314 mm (width x height x depth)
Weight	approx. 4.5 kg (without plug-in power supply)
Housing type of protection	IP 30
Electrical protective class	III
Test certificates	CE
Temperature allowed	Operation: +10 °C to + 35 °C (41 °F to 95 °F) Storage: -25 °C to + 65 °C (-13 °F to 268 °F)
Humidity allowed	≤ 80 %
Climatic class	2 / VDI/VDE 3540
Power pack	Type: EDACPOWER EA1036R Input: 100 - 240 V ~ / 50 - 60 Hz / 1 A Output: 12 V = / 3 A
EMC	EC directive 2004/108/EC EN 61326-1 – Interference emission: Class B – Interference immunity: IEC 61000-4-3 – Tolerance extension: 0.008 E FCC Class A
Meter safety	EC directive 2006/95/EC EN 61010-1

1.15 AVAILABLES LANGUAGES:

- German
- English
- Français
- Español
- Italiano
- Bulgarian/Български
- Česko
- Simplified Chinese/ 中文
- Traditional Chinese/ 繁體中文
- Greek/Ελληνικά
- Indonesian/Indonesia
- Japanese/ 日本語
- Magyar
- Malay/Melayu
- Norsk
- Polski
- Portugése
- Russian/Русский
- Slovenščina
- Thai/ ภาษาไทย
- Turkish/Turkce
- Dansk
- Română
- Nederlands

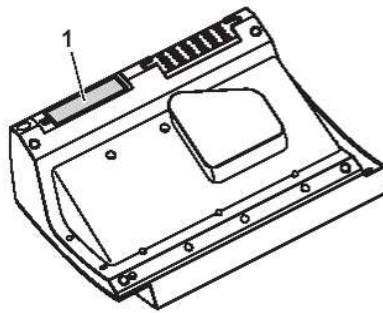
6 COMMISSIONING

For the first starting up, perform the following activities:

- Insert the buffer batteries
- Connect the power supply
- Switch on the photometer
- Set the language
- Set the date and time

1.16 BUFFER BATTERIES INSTALLATION

The buffer batteries supply the integrated clock while the photometer is switched off. Four alkaline manganese batteries (type AA or Mignon) separately included in the scope of delivery are used as the buffer batteries. Insert the batteries as follows:

1. Turn the photometer upside down and place it on a soft surface.
2. Open the lid of the battery compartment (1).
3. Insert the four batteries in the battery compartment. Make sure that the poles of the batteries are in the correct position.
4. The \pm signs on the batteries must correspond to the \pm signs in the battery compartment.
5. Close the lid of the battery compartment.

Batteries time life: The power consumption of the clock is very low. The lifetime of high quality batteries is at least 5 years.

1.17 CONNEXION OF POWER SUPPLY

The power is supplied via the enclosed plug-in power pack. The power pack supplies the photometer with low voltage (12 VDC).

CAUTION

The line voltage of the usage location must fulfil the specifications stated on the power pack. Always use the supplied 12 V original power pack only. Before plugging in the power cable check whether it is undamaged. If the power cable is damaged, the instrument must not be operated.

- 1) Connect the miniplug of the power pack to the socket (1) of the spectrophotometer.
- 2) Connect the power pack to an easily accessible power socket. The display illumination switches on and then off again.

1.18 CHOICE OF THE LANGUAGE

During the initial commissioning, the spectrophotometer automatically guides the user through the setting of the meter language, date and time after switching on (see following sections).

- Press key. <ON/OFF> the spectrophotometer is switched on.

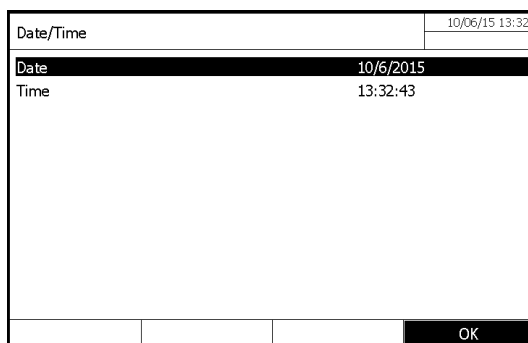
1. With «▲» and «▼» keys, select the language.
2. Confirm the choice with <START-ENTER> key.
The language has been set. The currently selected language is marked by a check.


NOTE: After the initial commissioning, the operator can change the language at any time. From the main menu <HOME> enter in the menu “General setup”, then “Language” menu.

1.19 DATE AND TIME SETTING UP

During the initial commissioning, the instrument automatically guides the operator to the setting of the time and date after the setting of the language.

The menu « Date / Time » is opened.

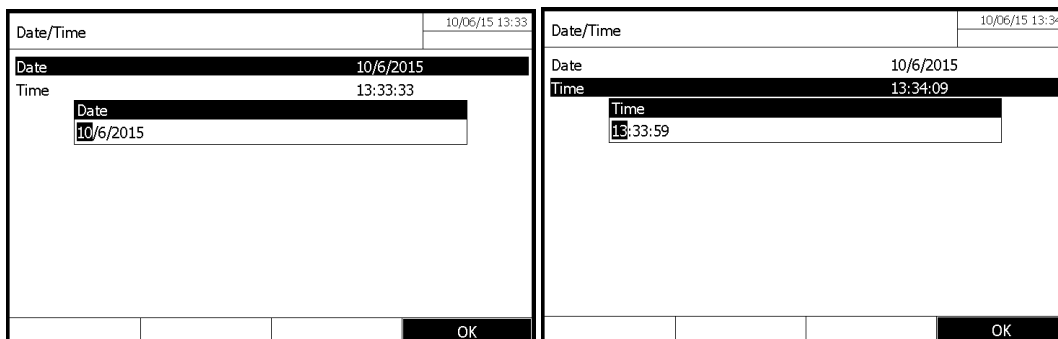1. With «▲» and «▼» keys, select one option from the menu. Valid this option with <START-ENTER> key.
2. The input field for the current date pops up.
3. Enter the current date with <0...9> keys and confirm.
4. The input field closes. The date is accepted.
5. Select « Time » and confirm.
6. The input field for the current hour pops up.

UviLine Series

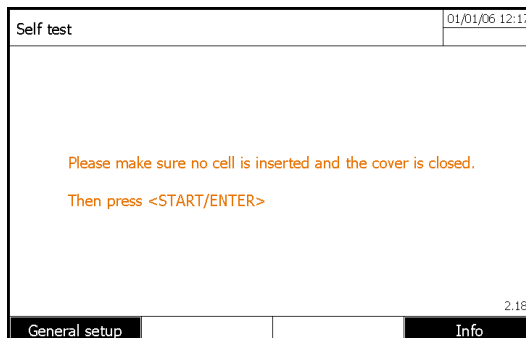
7. Enter the current hour with <0...9> keys and confirm.
The input field closes. The time is accepted.

NOTE: After the initial commissioning, the operator can change the date and time at any time. From the main menu <HOME> enter in the menu "General setup", then "Date/time" menu.


1.19.1 SELFT-TEST

IMPORTANT

During self-test, the cuvette compartment must be empty (without anymore cuvette inside the cuvette well) and the cover of the cuvette compartment must be closed.

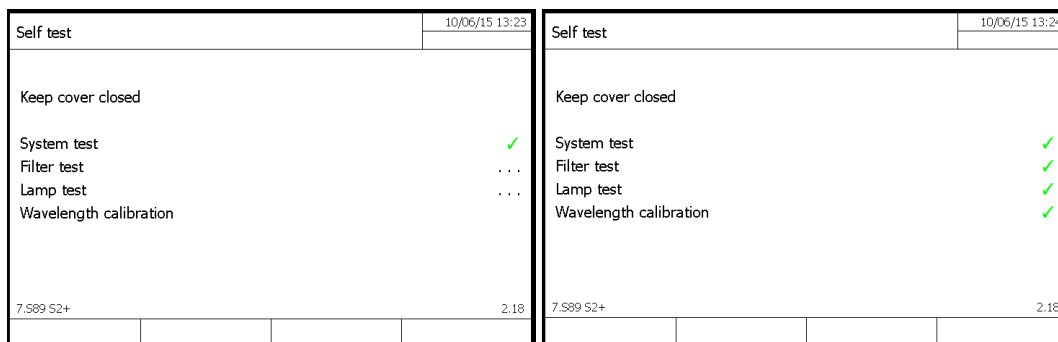- Press <START ENTER> key to run the self-test.

During self-test, different parts of the spectrophotometer are checked. If the test is OK, the symbol ✓ is displayed on the corresponding line.

This self-test includes:

- memory, processor, interfaces, filter and lamp tests.
- checking of the wavelength calibration.

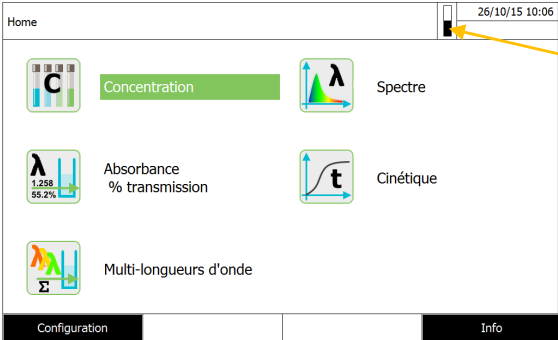
It is possible to see and print the self-test results by pressing F4 key [Info]

1.19.2 WARM-UP TIME

After switching on the photometer requires a warm-up time of 15 minutes. Reproducibility of measurement data is restricted during the warm-up time.

Therefore, do not measure during the warm-up time.


During the warm-up time, a progress bar appears on the display next to the date. The progress bar disappears as soon as the warm-up time is over.

Progress bar during warm-up time

7 OPERATION

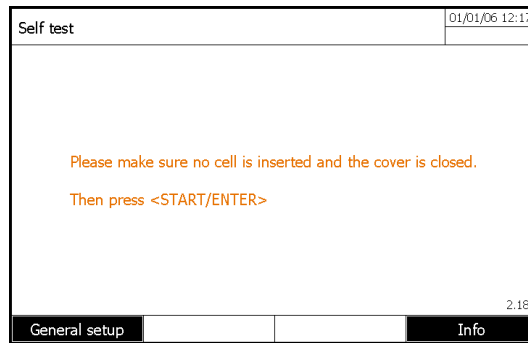1.20 SWITCHING ON OR OFF THE PHOTOMETER

Switch the photometer on with **<ON/OFF>**.

- The Self test dialog (if the user management is not active) or the Login dialog (if the user management is active)
- Enter user name and password or register as a guest then press <START-ENTER>

Switching off: To switch off the instrument, press the **<ON/OFF>** key.

1.21 GENERAL OPERATING PRINCIPLES

1.21.1 NAVIGATING

After the end of the self-test, the screen displays the main menu. 5 modes are available:

Press «▲» and «▼» keys, to select the wanted mode

- Concentration
- Absorbance / % Transmission
- Special / Multi wavelengths
- Spectrum
- Kinetics

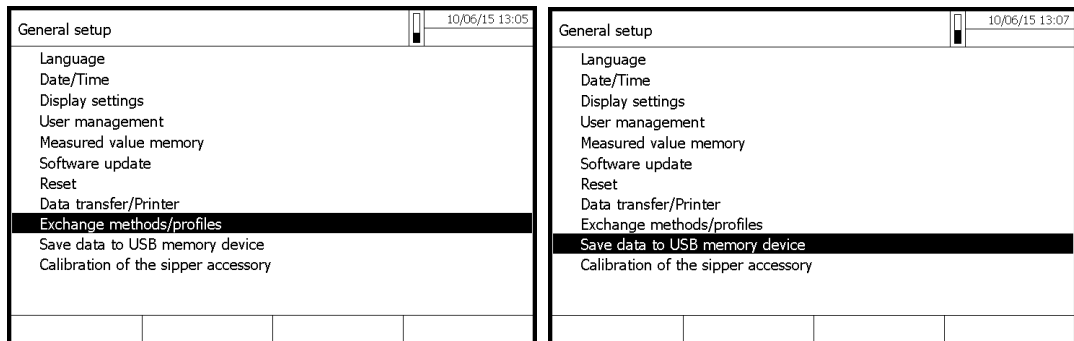
Press <F2> key from the main menu to access in general setup options

1.21.2 CUSTOMER METHODS AND PROFILE IMPORTATION/EXPORTATION

This function gives you possibility to export and import methods from an instrument to another one. Therefore, you can program all your methods in one instrument and duplicate them on all other same instrument easily with an USB stick.


Configuration menu:


You have the choice to either export all data of your instrument or to select on a list which emthods or profile you want to

1.21.3 PDF FILE PRINTING

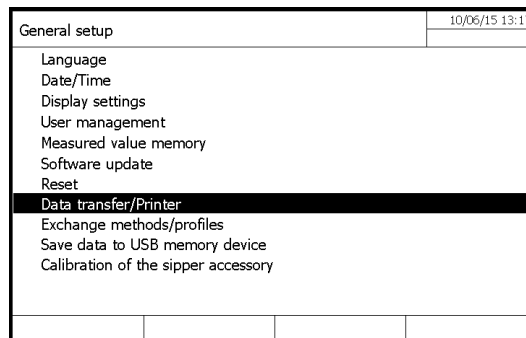This new function gives you possibility to print on a virtual printer by generating a printed report on a PDF file. This is useful when a physical printer is not available or you want to store numerically printed result directly.

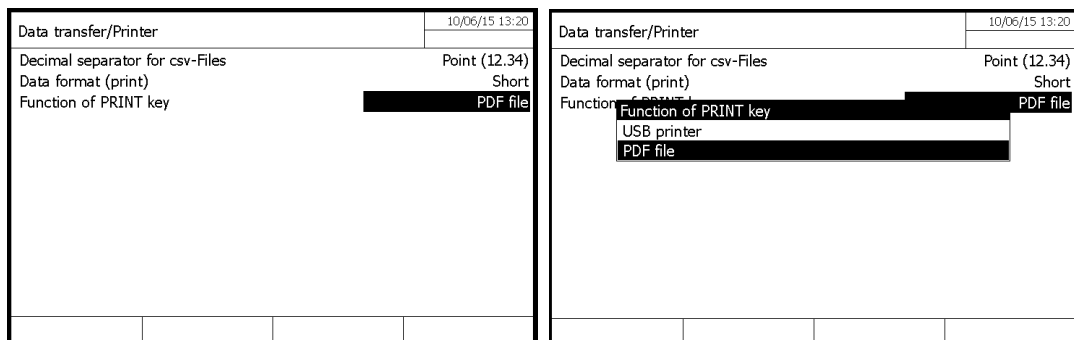General setup menu to configure printer parameters


With parameter Function of <PRINT> key, you can select if you want to use a real USB printer or the virtual printer generating PDF file printed report.

8 CONCENTRATION MODE



1.22 GENERAL POINT

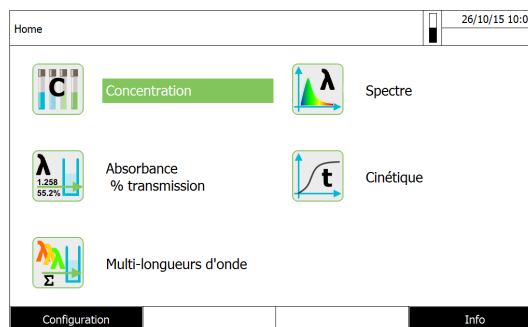The concentration displays results according to existing standard methods, or specific method defined by the user.

1.23 METHOD PROGRAMMING

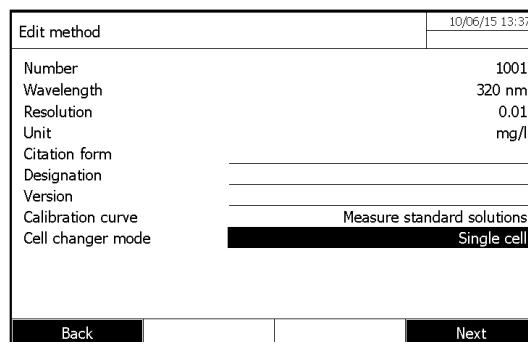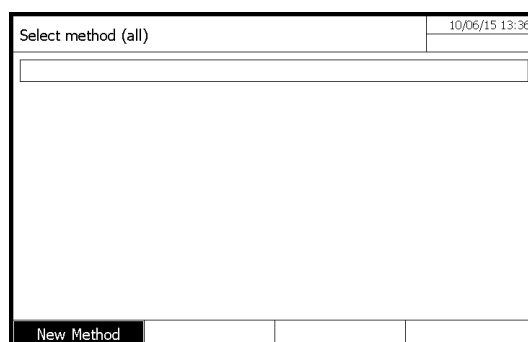In the "concentration" mode use F1 [New methods] for a new set up.

*** Number**

Automatic classification of methods (From 1001 to 1100).

*** Wavelength**

320 – 1100 nm for UviLine 8100C / 9100C / 9300

190 – 1100 nm for UviLine 9400C / 9600

*** Resolution**

From 0 to 3

It is the number of figures after comma for the result.

*** Unit**

i.e.: mg/L (10 characters maximum).

*** Citation form**

(18 characters maximum).

Enter the chemical formula of the compound to be measured. i.e.: NO₃

*** Designation**

(18 characters maximum).

Enter the compound name. i.e.: Nitrate

*** Version**

(18 characters maximum).

Enter the parameter concentration range. i.e.: 0.9 – 75.3 mg/L

*** Calibration curve**

There are several possibilities:

- Measure the standard solutions.
- Enter the equation of the calibration curve or the factor value.
- Simulate measurements while entering of the couples of Absorbance/Concentration values

*** Cell changer mode**

Usefull if the instrument is equiped with a cell changer

*** Standard ID**

Enter the identification number of the standards if this one exists.

*** Standard manufacturer**

Enter the standard manufacturer name if this one exists

*** Blank value**

YES or NO.: Measure done or not of the reagent blank

*** Standard conc 1**

Enter the concentration of standard 1.

UviLine Series

MEASURE OF STANDARD SOLUTIONS

Edit method		10/06/15 13:40
Standard ID	_____	
Standard manufacturer	_____	
Blank value	_____	No
Standard conc. 1	_____	10.00 mg/l
<div style="display: flex; justify-content: space-between;"> Back Add Next </div>		

Edit method		10/06/15 15:00
Standard ID	_____	
Standard manufacturer	_____	
Blank value	_____	No
Standard conc. 1	_____	10.00 mg/l
Standard conc. 2	_____	20.00 mg/l
Standard conc. 3	_____	30.00 mg/l
<div style="display: flex; justify-content: space-between;"> Back Add Delete Next </div>		

Remark

It is possible to add other standards (Maximum 10) by pressing F2 key [Add] and to enter their concentration. To delete a standard, select it with «▲» and «▼» keys and press F3 key [Delete]. When all standards are entered, press F4 key [Next] to continue.

ENTER VALUE PAIRS

Edit method		10/06/15 15:02
Number	1001	
Wavelength	320 nm	
Resolution	0.01	
Unit	mg/l	
Citation form	_____	
Designation	_____	
Version	_____	
Calibration curve	_____	Enter value pairs
Cell changer mode	Single cell	
<div style="display: flex; justify-content: space-between;"> Back Next </div>		

Edit method		10/06/15 15:03
Standard ID	_____	
Standard manufacturer	_____	
Blank value	_____	Yes
Standard conc. 1	_____	10.00 mg/l
Standard conc. 2	_____	20.00 mg/l
Standard conc. 3	_____	30.00 mg/l
<div style="display: flex; justify-content: space-between;"> Back Add Delete Next </div>		

ENTER FORMULA

Edit method		10/06/15 15:03
Number	1001	
Wavelength	320 nm	
Resolution	0.01	
Unit	mg/l	
Citation form	_____	
Designation	_____	
Version	_____	
Calibration curve	_____	Enter formula
Cell changer mode	Single cell	
<div style="display: flex; justify-content: space-between;"> Back Next </div>		

Edit method		10/06/15 15:03
$c = a_0 + a_1 \cdot A + a_2 \cdot A^2 + a_3 \cdot A^3 + a_4 \cdot A^4 + a_5 \cdot A^5$		
a0	_____	
a1	_____	
a2	_____	
a3	_____	
a4	_____	
a5	_____	
Lower limit of measuring range	_____	
Upper limit of measuring range	_____	
<div style="display: flex; justify-content: space-between;"> Back Next </div>		

Enter a function as: $C = a5.A^5 + a4.A^4 + a3.A^3 + a2.A^2 + a1.A + a0$

With **C** = Concentration
A = Absorbance
a0 to **a5** free factors (from 0.000 to 9999.000)

Example: Use of a K factor K (C = a1A)
 - a0 = 0
 - a1 = K (factor)
 - a2, a3, a4, a5 = Zero

Example: straight line regression (C = a1A + a0)
 - a0 = bias
 - a1 = slope
 - a2, a3, a4, a5 = Zero

Example: third degree curve (Nitrate test in sea water) : $C = 0.1493A^3 - 2.5154A^2 + 36.524A + 1.865$
 - a0 = 1.865
 - a1 = 36.524
 - a2 = - 2.5154
 - a3 = 0.1493
 - a4 & a5 = 0

Edit method	10/06/15 15:04
c = a0 + a1*A + a2*A ² + a3*A ³ + a4*A ⁴ + a5*A ⁵	
a0	_____
a1	_____
a2	_____
a3	_____
a4	_____
a5	_____
Lower limit of measuring range	0.00 mg/l
Upper limit of measuring range	75.00 mg/l
Back	Next

*** Lower and upper limit of measuring range**
 Choice between zero and the maximum value of measured parameter.

TIMER

If your methods needs timers, you can edit vaues by pressing <F4>, [NEXT]

Edit method	01/01/15 12:02
Timer 1	00:00:00
Timer 2	00:00:00
Timer 3	00:00:00
Timer 4	00:00:00
	0:25:00
Back	Complete

1.24 METHOD SAVING

The methods are automatically saved under the classification name from 1001 to 1100.

Remark: However, methods using standards will be saved only after standards measurement.

1.25 METHOD DELETING

- Select method to be deleted in the method List with «▲» and «▼» keys and press <START/ENTER> key. Then press F1 key [Setup].
- Select « Delete method » with «▲» and «▼» keys and press <START/ENTER> key.

Concentration		10/06/15 15:05	
Zero measurement required! Press <ZERO/BLANK>			
1001		0.00 - 100.00 mg/l	
Setup	Method list		

Concentration		10/06/15 15:06	
Dilution			
Sample blank value			
✓ Display absorbance			
New method			
Edit method			
Delete method			
Measurement data memory			

1.26 METHOD EXECUTION

1.26.1 WITH STANDARD SOLUTIONS

1.26.1.1 Measure

When standard (s) are entered, press F4 key [Next]. Then - press F3 key [Measurement] to start the measure.

Edit method		04/16/15 13:13	
Blank value			
Std. 1	10.00 mg/l		0.000
Std. 2	20.00 mg/l		
Std. 3	30.00 mg/l		
Back	Measurement	Next	

Blank value ()		04/16/15 13:17	
Zero measurement required! Press <ZERO/BLANK>			
Nitrate			

- Press on <ZERO/BLANC> key to perform the zero of the meter.

Blank value ()		04/16/15 13:18	
Start measurement with <START/ENTER>			
Nitrate			

Blank value ()		04/16/15 13:19	
Last measured absorbance 0.377			
Median 0.377 (1 Measurement(s))			
Nitrate			
Next meas.	Discard		Apply

- Install the blank (if necessary) inside the cuvette holder.
- Start the blank measurement (if this one was programmed) by pressing <START/ENTER> key.
- It is possible, in the event of error, to refuse the value and to measure again the solution. Press then on F2 key [Discard].
- Then remake measurement while pressing on <START/ENTER> key.

- It is also possible, if the solution is unstable, to measure it again several times. The software will make the average of the various measured values.

Remark: The number of times to measure again the solution is not limited.

Blank value ()	04/16/15 13:20
Last measured absorbance 0.379	
Median 0.377 (4 Measurement(s))	
Nitrate	
Next meas.	Discard Apply

Blank value ()	04/16/15 13:20
Apply Last measured absorbance Median	
Median 0.377 (4 Measurement(s))	
Nitrate	
Next meas.	Discard Apply

- Press, then F4 key [Apply].

- The operator can choose to preserve the last measured value by validating the option "Last measured absorbance" or to use the average value calculated by choosing and validating the option "Median".

Edit method	04/16/15 13:22	
Blank value	10.00 mg/l	0.377
Std. 1	20.00 mg/l	0.000
Std. 2	30.00 mg/l	
Std. 3		
Back Measurement Next		

Std. 1 (10.00 mg/l)	04/16/15 13:23
Start measurement with <START/ENTER>	
Nitrate	

- Install the first standard to be measure. Then press F3 key [Measurement].

- Start measurement of the first standard by pressing <START ENTER> key.

- As for the blank, the operator can reject measurement (F2 key [Discard]) or make several times the measurement of the same standard (F1 key [Next meas.]) and preserve the average value.

- Make, in the same way, the measurement of the other standards.

Std. 1 (10.00 mg/l)	04/16/15 13:23
Last measured absorbance 0.884	
Median 0.884 (1 Measurement(s))	
Nitrate	
Next meas.	Discard Apply

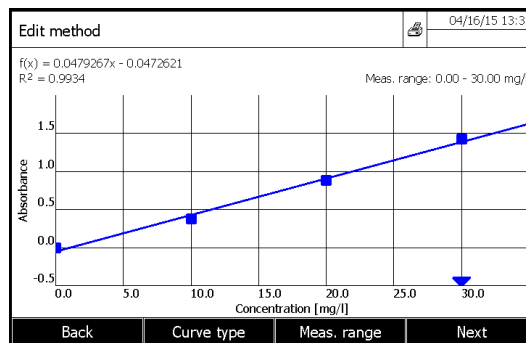
Edit method	04/16/15 13:24	
Blank value	10.00 mg/l	0.377
Std. 1	20.00 mg/l	0.884
Std. 2	30.00 mg/l	1.700
Std. 3		2.753
Back Measurement Next		

- Press F4 key [Next.] to proceed

NOTE:

If the programming of standards concentration is not increasing or decreasing in a monotonous way or if the measurement of the blank and standards absorbances is not, also increasing or decreasing in a monotonous way (see example below), the spectrophotometer will display an error prompt (see below)

1.26.1.2 Calibration curve displaying


F1 key [Back]

- It allows to the user, in the event of error, to return back to the analysis programming and to modify the acquiring way of the values to make the curve (standards measurement, use of the curve equation or entry of the Concentration/absorbance couples). That also enables him to change the standards concentration.

F2 key [Curve type]

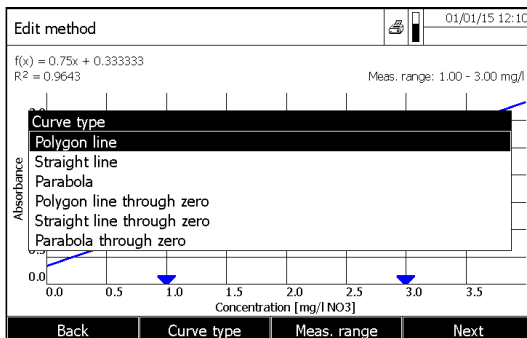It allows changing the curve calculation mode.

Straight line

The straight line calibration curve is consisted of line segments connecting the calibration points between them.

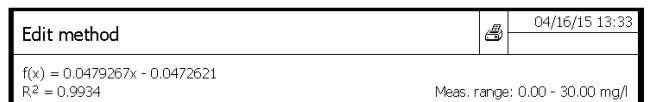
Linear regression

Curve whose equation is form $ax + B$.

Quadratic regression

Curve which equation is: $ax^2 + bx + c$.

The equation of the calculated curve is displayed on the top with the coefficient of R^2 determination

F3 key [Meas.range]

It allows to the operator to reduce the parameter measuring range.

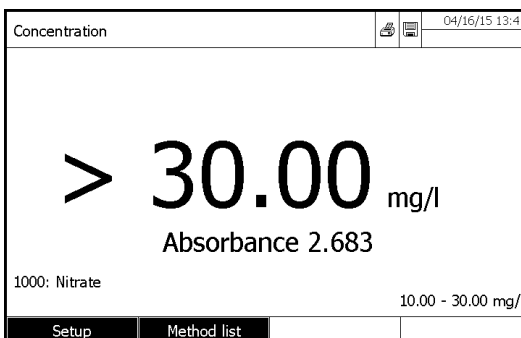
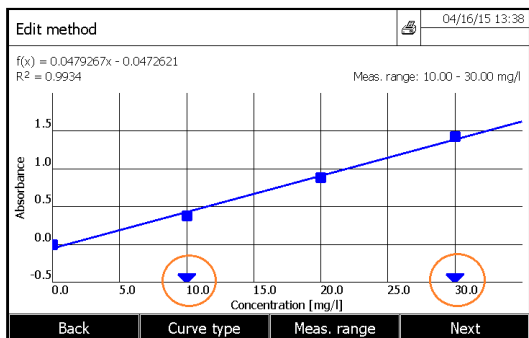
Lower limit
10.00 mg/l

Lower limit entered by the operator

Upper limit
30.00 mg/l

Upper limit entered by the operator



These limits are materialized by an arrow «▼» on the concentration axis.

Any value under or above limits fixed by the operator will be declare as > or < to the limits, but absorbances will be displayed (see example above).

1.26.1.3 Point deleting

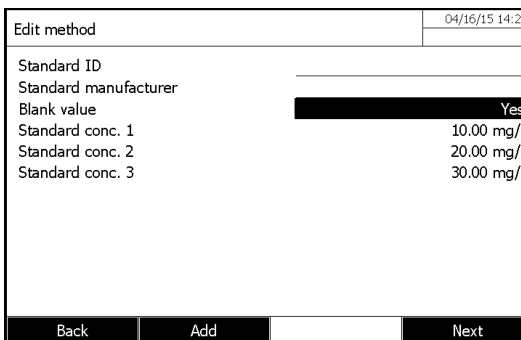
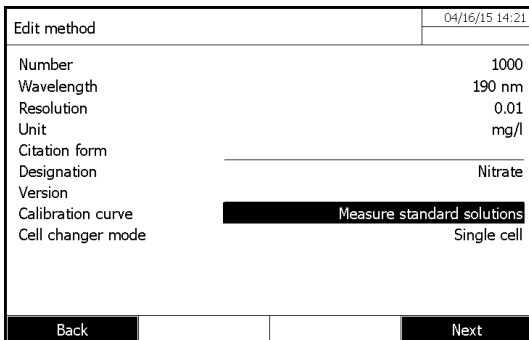
To delete a point press F1 key [Back], then Press F4 key [Next].

-- Select the point to delete with «▲» and «▼» keys and delete the point by pressing F3 key [Delete]. Then press on F4 key [Next].

During this stage, it is always possible to measure a standard once again.

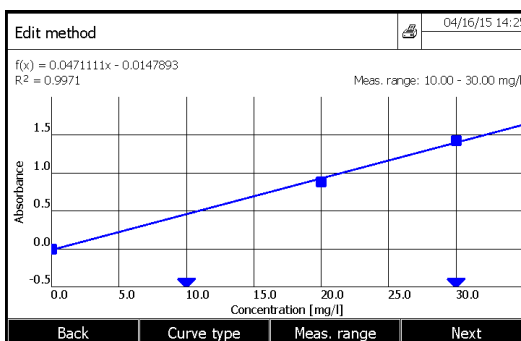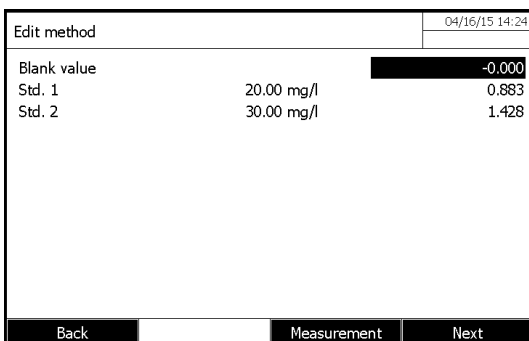- Select the standard to measure again.
- Press F3 key [Measurement] to start the standard measurement.
- Press F4 key [Next] to display the curve.


- To set back the removed standard, press F4 key [Back].
- Add the standard.
- Enter its concentration.
- Measure the concentration of this standard once again, or enter its absorbance following the curve building mode selected.

UviLine Series


1.26.1.4 Calibration curve printing

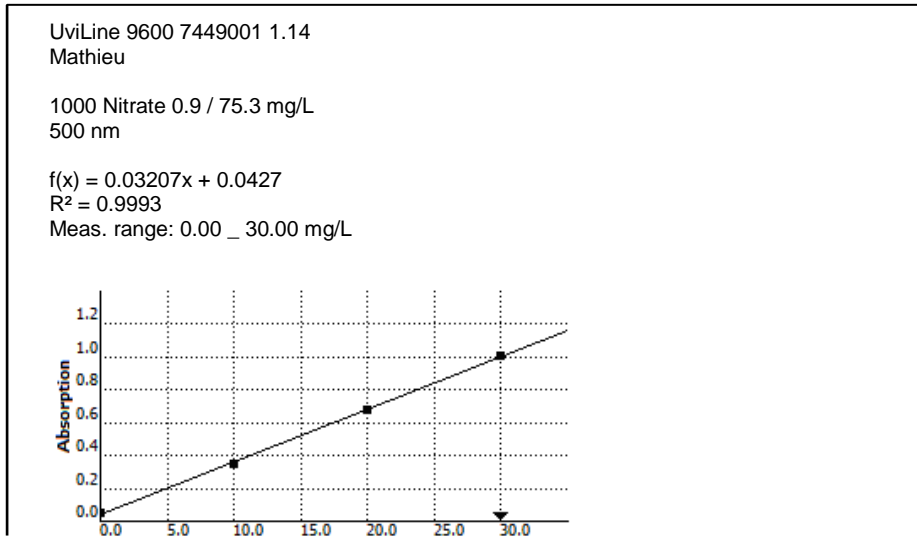When the curve is displayed, press the <PRINT> key.


1.26.1.5 Sample measurement

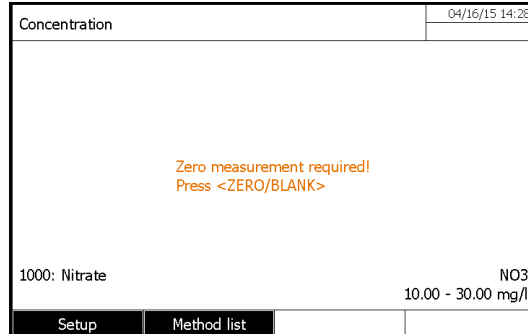
- Press <ZERO BLANK> key to start the blank measurement. It performs the zero

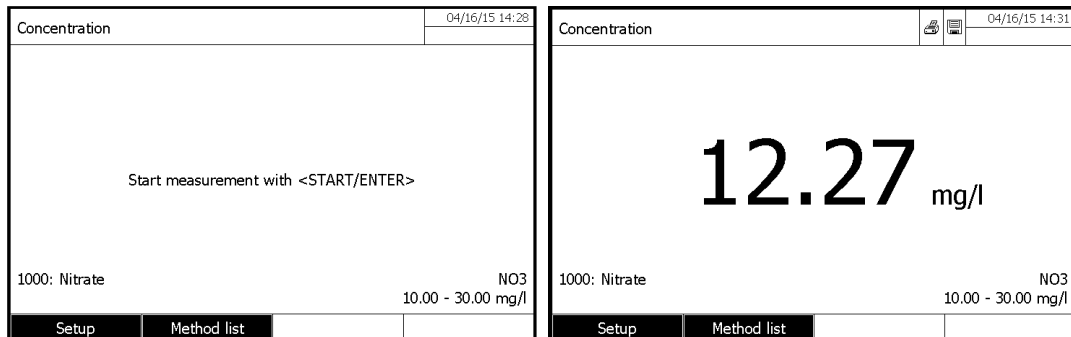- Press < START ENTER> key to start sample measurement.

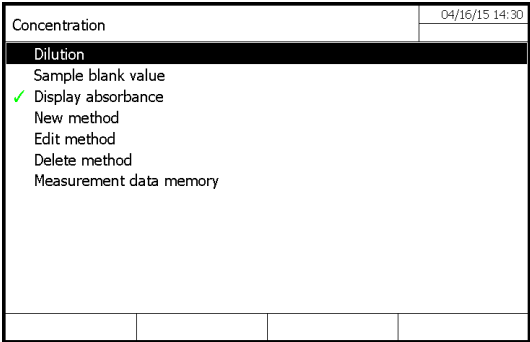
If you want that the samples absorbance is displayed at the same time as the concentration, he has to select it in the [Setup] menu and unselect it. The symbol ✓ means that the absorbance will be displayed at the same time as their concentration.


1.26.1.6 Results printing

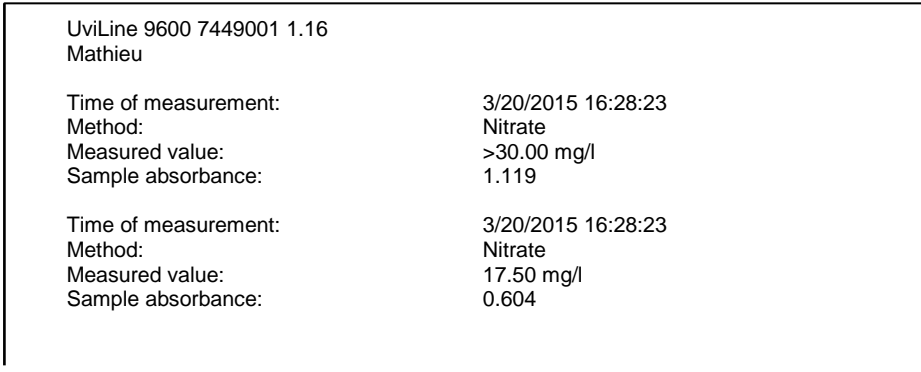- Press <PRINT> key to send data on printer.


1.26.1.7 Sample dilution

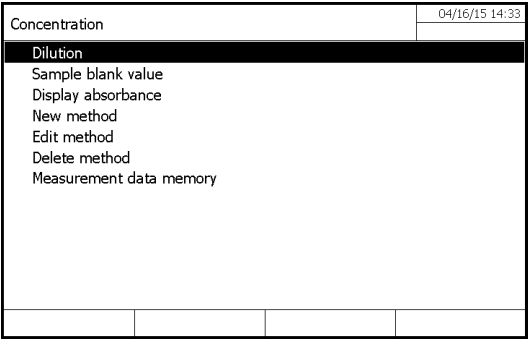It is possible to program a dilution factor (From 0 to 999) that will be taken in account during sample measurement. Before sample measurement. Press F1 key [Setup]. Then press on <START ENTER> key.

Enter the dilution factor.
1 + 0 = No dilution.
1 + 1 = Dilution by 2.

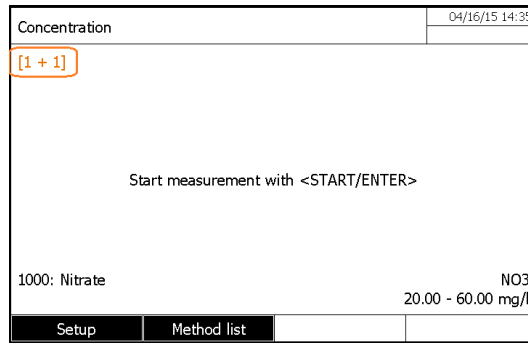
When the dilution facto is entered, confirm with <START ENTER> key. The dilution factor (i.e.: [1 + 1]) is displayed on the top left part of the screen.

NOTE: This dilution factor will be maintained with the value entered by the operator for all measurements to come. At the time of the exit of the mode of analysis or analysis in progress it will be given to 0 (no dilution).


1.26.1.8 Sample blank value

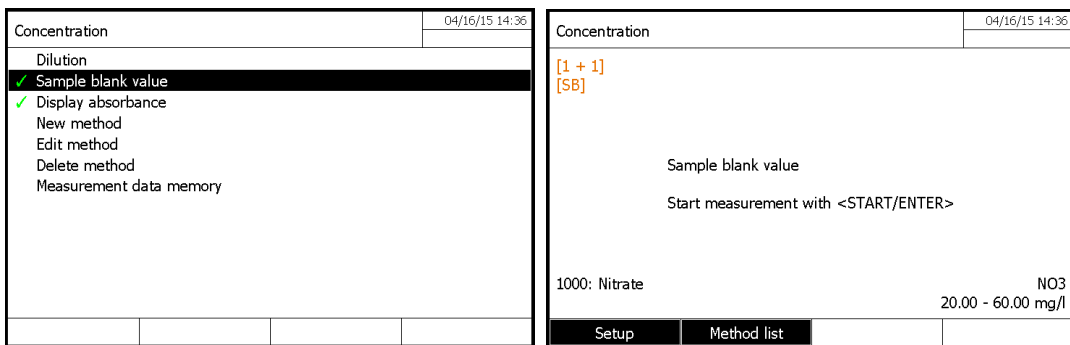It is possible to subtract a sample blank for each measured sample. Press F1 key [*Setup*] after zero measurement. It displays:


Remark: The symbol ✓ at the end of « Sample blank value » means that the option is selected.

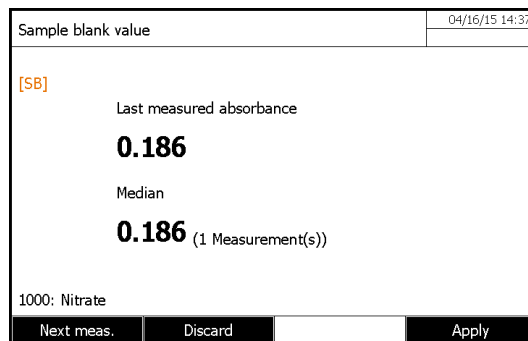The symbol [SB] appears at top left corner of the screen and stays here and will remain all the time that this option is selected or that the operator did not leave the analysis in progress or the mode of concentration measurement. Insert the first sample blank and run the measurement by pressing <START ENTER> key.

As for the blank, the operator can reject measurement (F2 key [*Discard*]) or make several times the measurement of the same sample blank (F1 Key [*Next meas.*]) and preserve the average value.

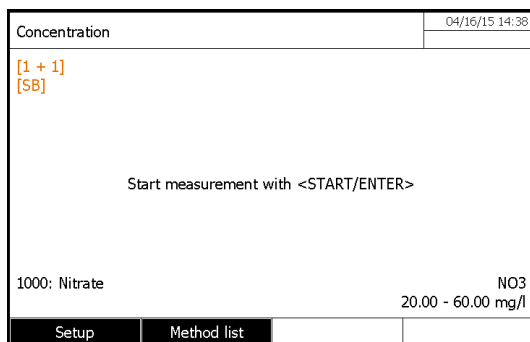
- Press F4 key [*Apply*] to run the sample measurement:
- Insert the sample corresponding to the measured sample blank and start the measurement by pressing <START ENTER> key.

It displays the sample absorbance read against the zero (i.e.: 0,562) and the sample concentration calculated as following:


Sample concentration = sample absorbance read against the zero minus sample blank absorbance read against the zero multiply by the factor or reported to the calibration curve.

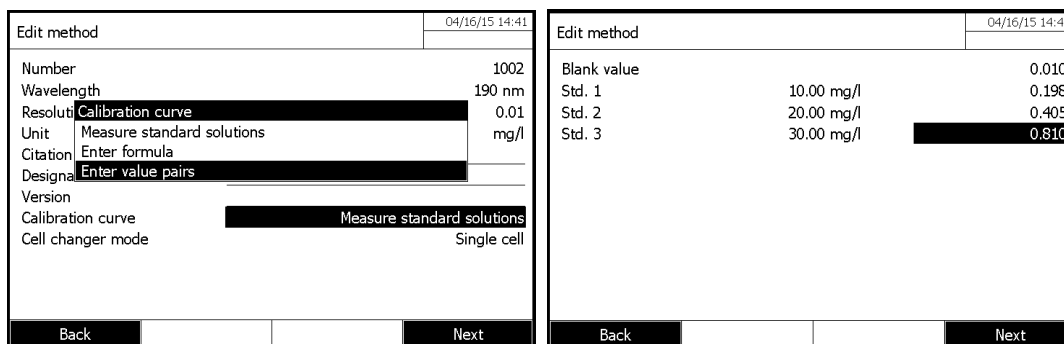1.26.2 WITH VALUE PAIRS

Enter, with keyboard, the absorbance values of the blank (if necessary) and the one of each standard and press F4 key [Next]. It displays then the calibration curve.

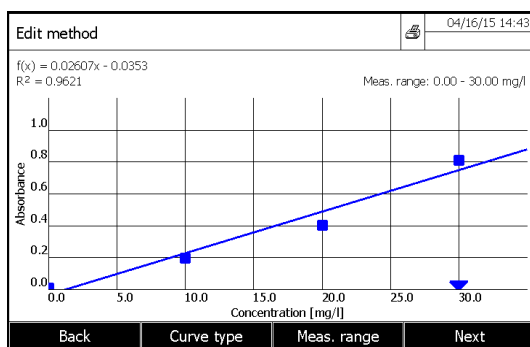From this point, the continuation proceeds is the same as previous chapter "standards.Solutions"

1.26.3 WITH A FORMULA

Edit method		04/16/15 14:44
Number		1002
Wavelength		190 nm
Resolution	Calibration curve	0.01
Unit	Measure standard solutions	mg/l
Citation	Enter formula	
Designation	Enter value pairs	
Version		
Calibration curve	Enter value pairs	
Cell changer mode		Single cell

Edit method		04/16/15 14:46
c = a0 + a1·A + a2·A ² + a3·A ³ + a4·A ⁴ + a5·A ⁵		
a0		
a1		
a2		
a3		
a4		
a5		
Lower limit of measuring range		
Upper limit of measuring range		

When the formula is entered, press F4 key [Next], then proceed as previous chapter "standards.Solutions"

Concentration		04/16/15 14:48
Zero measurement required! Press <ZERO/BLANK>		
1002		0.00 - 100.00 mg/l

1.27 METHOD EDITING

- Enter in « Concentration » mode.
- Select a method with «▲» and «▼» keys, then confirm with <START ENTER> key.
- Press F1 key [Setup].

Concentration		04/16/15 14:48
Zero measurement required! Press <ZERO/BLANK>		
1002		0.00 - 100.00 mg/l

Concentration		04/16/15 15:10
Dilution		
Sample blank value		
✓ Display absorbance		
New method		
Edit method		
Delete method		
Measurement data memory		

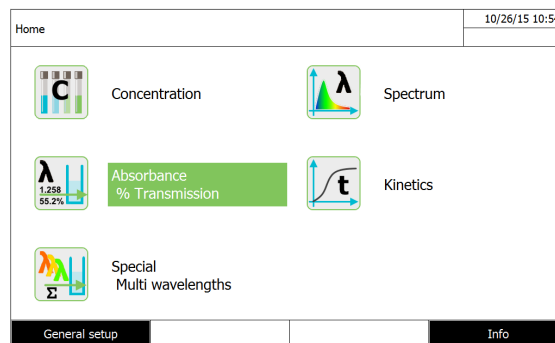
- Select « Edit method » with «▲» and «▼» keys, then confirm with <START ENTER> key.

9 ABSORPTION / TRANSMISSION MODE

1.28 GENERAL POINTS


The absorbance or transmittance respectively is measured without the use of any methods. All settings are configured during measurement.

The absorbance or transmittance can alternatively be measured against the air or against a reference solution determined by the operator. The settings of the last measurement are active.

1.29 MEASUREMENT

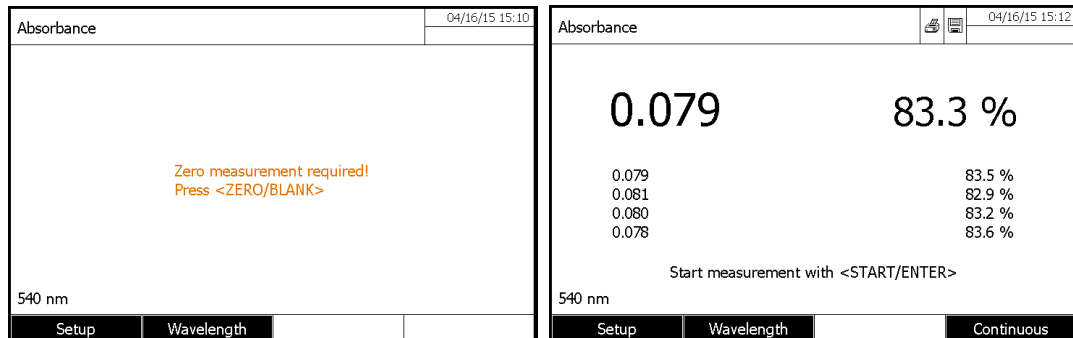- Press F2 key [*Wavelength*] to enter a new wavelength.
- Press <ZERO BLANK> key to start the zero measurement.
- Press <START ENTER> key to start the sample measurement (Absorbance and transmittance are displayed).
- Press <START ENTER> key to start a new measurement.


When there were several measurements, the four last are displayed at the same time as measurement in progress

- Press F4 key [*Continuous*] to measure the same sample several time again.
- Press F4 key [*Stop*] to stop the continuous measurement.

1.30 RESULTS PRINTING

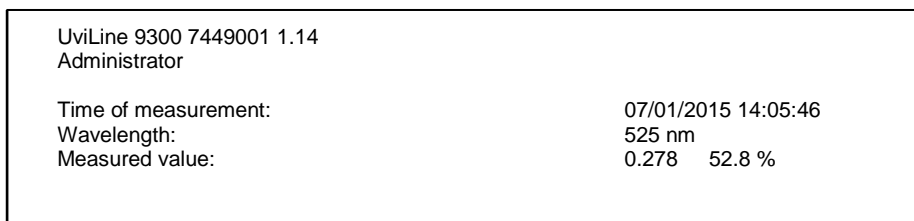- Press <PRINT> key to print the results.


10 SPECIAL/ MULTI-WAVELENGTHS

1.31 BASIC INFORMATION

In the Special / Multi wavelengths mode of the UviLine Series, you can carry out measurements with special methods and functions. You can use the following functions for these methods:

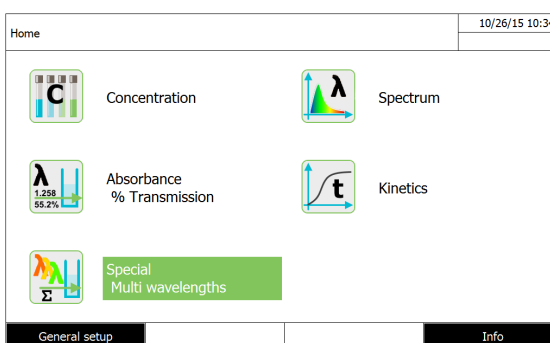- Measurements at different wavelengths
- Use of procedure variables. Procedure variables provide a value that has to be entered prior to each measurement on the photometer (e.g. volume, pH value or temperature)
- Check whether a value meets a condition. With a condition you can check a value for validity (e.g. absorbance value, procedure variable or the result of a formula).
- Formula editor for the convenient programming of any user-defined methods

1.32 METHOD PROGRAMMING

Remark

For multi wavelength methods, you can use the method numbers 2001 to 2499. All special methods can also be selected in the method list of the concentration mode.

The creation of a user-defined method is done in the following steps:

- **Enter the general method data**
Method number, method name, version, unit etc.
- **Enter the wavelengths for absorbance measurements (Ax nm)**
Minimum 1, maximum 10
- **Define the procedure variables (KX) (optional)**
Procedure variables are used to take into account any influence quantities that cannot be measured by the photometer.
The values for these procedure variables have to be entered for all measurements with the method, e.g. the temperature or pH value.
- **Enter the formula to calculate the measurement result**
Enter the formula with which you want to calculate the measurement result in the formula editor.
- **Enter an additional condition (optional)**
Conditions are used to check the measurement result for validity.
The condition is entered with the formula editor.

To create a new method:

- Press F1 key [*New method*].
- Enter here general data of the method. The next number of method available is already registered as next number of analysis to be used.

To fill the fields with entry, it is possible to proceed as follows:

- Fill the entire empty fields with entry ones after the others.
- Press on the key F2 [*Method list*] to select a method already recorded
- Press on the F3 key [*Delete*] to remove the method completely.
- Press on the F4 key [*Next*] to validate all the entries and to pass in the following page.

Edit method		01/01/15 12:09
Number	2002	
Name		
Version		
Citation form		
Unit		
Resolution	0.01	
Lower limit of measuring range		
Upper limit of measuring range		
Cell changer mode	Single cell	
Method list		Delete Next

Edit method		01/01/15 12:30
Number	2055	
Name	Protein	
Version	1.1	
Citation form	Prot	
Unit	mg/mL	
Resolution	0.01	
Lower limit of measuring range	1.00 mg/mL	
Upper limit of measuring range	100.00 mg/mL	
Cell changer mode	Single cell	
Method list		Delete Next

- Press on F2 key [Add] to add another wavelength.
- Press F3 key [Effacer] to delete the last wavelength.
- Press on the F4 key [Next] to validate all the entries and to pass in the following page.
- Add variable if necessary (for example 'sample temperature')
- Press on the F4 key [Next] to validate all the entries and to pass in the following page.

Wavelength		01/01/15 12:15
Wavelength 1	280 nm	
Wavelength 2	260 nm	
Back		Add Delete Next

Procedure variables		01/01/15 12:32
K1	Temp	
K2	Pressure	
Back		Add Delete Next

- Start writing the formula by using operators (F2) and variables (F3)
- Press on the F4 key [Next] to validate all the entries and to pass in the following page.

Formula entry		01/01/15 12:36
<p>Use the <Operators> softkey to select an operation, function or constant (e.g.: +, -, *, tan, log, e, Pi).</p> <p>Use the <Variables> softkey to select an absorbance at a certain wavelength or a procedure variable.</p> <p>Enter numerals via the keyboard.</p> <p>You can erase the last entry with <<>.</p>		
Back		Operators Variables Next

Formula entry		01/01/15 12:38
$R = (A_{230 \text{ nm}} \cdot 0.183 - 0.075 \cdot A_{260 \text{ nm}}) \cdot K_1 / K_2$		
Back		Operators Variables Next

- You can define a condition formula wish validate or not the measure (for example: limit absorbance value)

Condition		01/01/15 12:39
<p>Here you can enter a formula for a condition. The measured value is only valid if this condition is met.</p>		
Back		Operators Variables Next

Condition		01/01/15 12:44
$A_{230 \text{ nm}} < 3.1$		
Back		Operators Variables Next

- Press on the F4 key [Next] to save and start the method.

1.33 METHOD DELETING

- Select the method to be deleted in the multi-wavelength mode with «▲» and «▼» keys and confirm with <START ENTER> key.
- Press F1 key [Setup].
- Select « Edit method » with «▲» and «▼» keys and confirm with <START ENTER> key.

Edit method		01/01/15 12:46
Number	2000	
Name	test	
Version	1	
Citation form		
Unit		
Resolution	0.01	
Lower limit of measuring range		
Upper limit of measuring range		
Cell changer mode	Single cell	
Method list		Delete Next

- Press F3 key [Delete].

1.34 METHOD SELECTION

Select method (all)		01/01/15 12:50
<input type="text"/>		
2055	Protein	Prot 1.00 - 100.00 mg/mL
		Last used

The method list is displayed.

Methods are ordered by method number.

Method selection:

- Select a method with «▲» and «▼» keys. The active selection is displayed in reverse video.

LIMITATION OF METHODS LIST

If the list is very long, the operator can narrow down the method list and thus make the search easier as follows:

- Press F4 key [Last used], it is possible to restrict the method list to the ten methods last used.
 - With the search function, the operator can search certain character strings in the list. The search takes place as a full-text search of the entire list contents.
- The operator can search for a method number or certain citation form.

Search for a character string:

- Enter the character string to be searched for in the search window with <A...9>.

The list appearing below shows all hits containing the character string. The hit list is updated with each character that is entered.

Select method (all)		01/01/15 12:51
<input type="text" value="Pro"/>		
6 m n o M N O		
2055	Protein	Prot 1.00 - 100.00 mg/mL
		Last used

1.35 METHOD EXECUTION

- Select a method with «▲» and «▼» keys and confirm with <START-ENTER> key in the multi-wavelength mode.
- Enter de variables values by using F4
- Press <ZERO BLANK> key

Special / Multi wavelengths		01/01/15 12:53		Procedure variables		01/01/15 12:55	
Zero measurement required! Press <ZERO/BLANK>				K1: Temp 22			
				K2: Pressure 1.01			
2055: Protein		1.00 - 100.00 mg/mL		Prot			
Setup	Method list	Transmission	Variables	Back			Apply

- Zero is performed at the programmed wavelengths.
- Press <START ENTER> key to start the sample measurement.

Special / Multi wavelengths		01/01/15 12:56		Special / Multi wavelengths		01/01/15 12:58	
Start measurement with <START/ENTER>				A(230 nm) = 0.919 A(260 nm) = 0.958			
				2.10 mg/mL			
2055: Protein		1.00 - 100.00 mg/mL		Prot			
Setup	Method list	Transmission	Variables	Setup	Method list	Transmission	Variables

It displays at the same time the absorbance values (or transmittance values) of the sample read at the different wavelengths.

- Press on F3 key [*Transmission*] to have the transmission values (%).
- Press on <START ENTER> key to run a new measurement.
- If the condition is not met, folowing message is displayed

Special / Multi wavelengths		01/01/15 12:57	
A(230 nm) = >3.300		A(260 nm) = >3.300	
Condition not met - - - - mg/mL			
2055: Protein		Prot 1.00 - 100.00 mg/mL	
Setup	Method list	Transmission	Variables

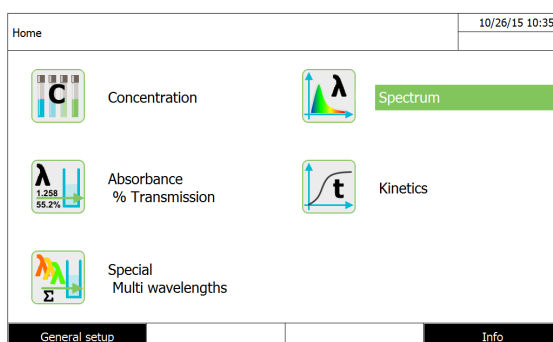
11 SPECTRUM MODE

1.36 GENERAL POINTS

The Spectrum mode makes it possible to measure and record the Absorption and Transmission values according to the wavelength. The wavelength range can be freely selected inside the spectrophotometer measuring range.


Baseline

A baseline has to be recorded before a spectrum is recorded. The baseline has to cover at least the wavelength range of the spectrum to be recorded. Once the baseline is measured, it remains stored in the photometer until a new baseline is recorded, the *Spectrum* mode is exited or the photometer is switched off.

1.37 METHOD PROGRAMMING

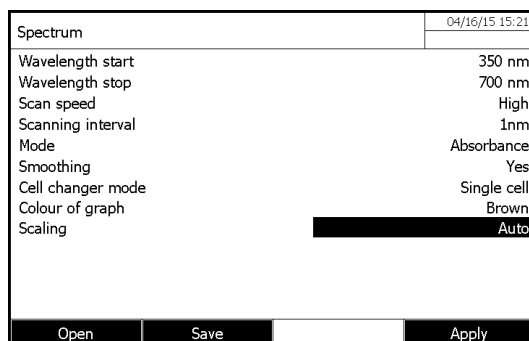In the "spectrum" mode, press on F1 key [Setup].

* Start wavelength

From 320 nm for the UviLine 8100C / 9400C / 9300
From 190 nm for the UviLine 9400C / 9600

* Stop wavelength

Up to 1100 nm for all UviLine

* Scan speed

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* Scanning interval

1, 2, 5, or 10nm interval point resolution.

* Mode

Absorbance or Transmission

* Smoothing

YES or NO

Cell changer mode

YES or NO

* Colour of the graph

Brown; Blue; Orange; Green; Red; Purple

* Scaling

Auto or Manual

1.38 METHOD SAVING

- Enter in the « Spectrum » mode.
- Press on F1 key [Setup] to enter inside the programming parameters sub-menu.
- Program parameters.
- Press on F2 key [Save].
- Press on F1 key [Location].
- Select the place where the method will store (Internal folder DataB or USB memory)
- Give a name to the method and confirm with <START/ENTER> key.

1.39 METHOD DELETING

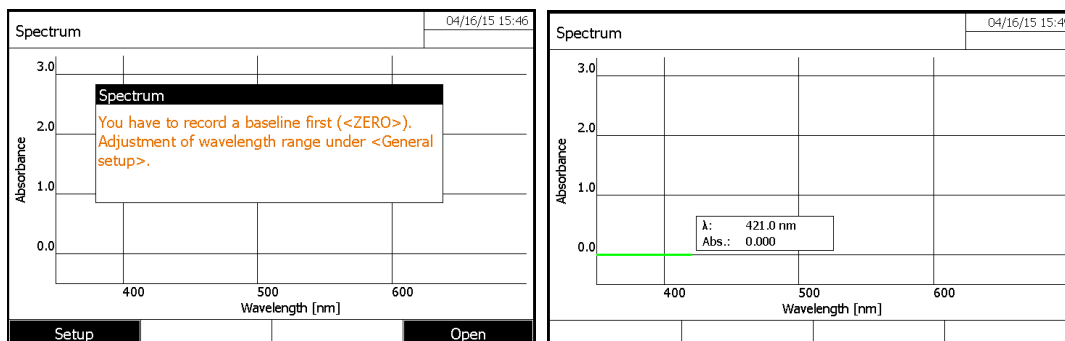
- Enter in the « Spectrum » mode.
- Press on F1 key [Setup].
- Press on F1 key [Open]. It displays:
- Select the method to be deleted with «▲» and «▼» keys.
- Press on F2 key [Delete].

1.40 METHOD SELECTION


- Enter in the « Spectrum » mode.
- Press on F1 key [Setup] to enter inside the programming parameters sub-menu.
- Press on F1 key [Open].
- Press on F1 key [Location] if the method is stored inside a USB memory.
- Select the method and confirm with <START/ENTER> key.

1.41 MEASUREMENT EXECUTION

- Enter in the « Spectrum » mode.
 - Start the baseline measurement by pressing <ZERO BLANK> key. The spectrophotometer memorizes the baseline.
- Wait until the end of measurement.

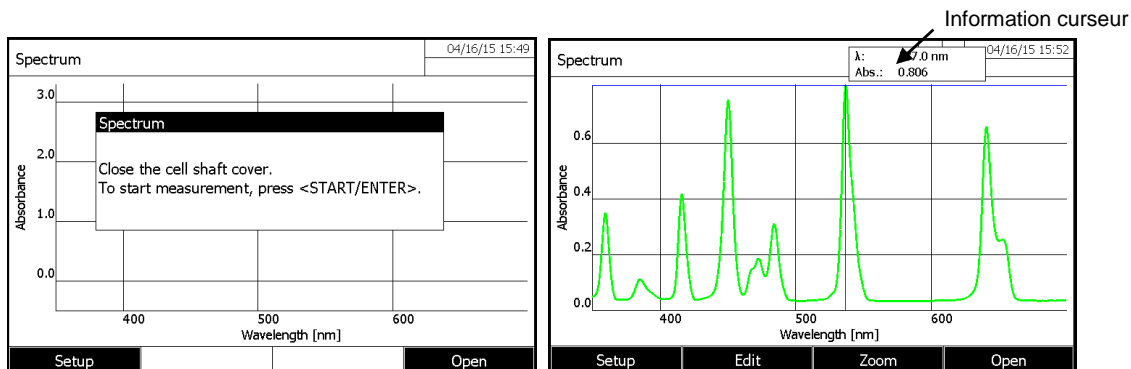- Insert the sample inside the cuvette holder, close the cover the ress <START ENTER> key to start the measurement.

The cursor appears at the absolute maximum of the spectrum.

The operator has the following possibilities:

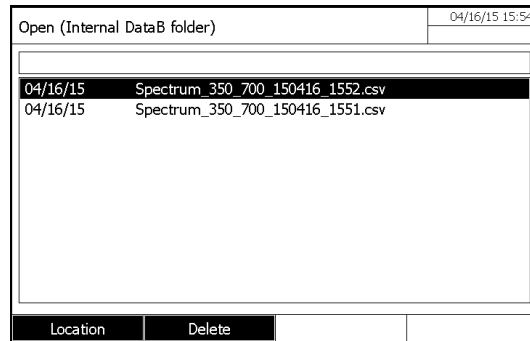
- Immediately realized some calculations on the spectrum (see following paragraph).

- To print, with <PRINT> key, the spectrum on a connected printer as a graphic.
- To record, with <STORE> key, the spectrum as a "*.csv" file. As the memory location, it is possible to select the spectrophotometer (*Internal DataB folder*) or a connected USB memory at the USB-A connection (*USB memory*).

1.42 SPECTRA TREATMENT


1.42.1 DOWNLOADING OF STORED SPECTRUM

Enter in the « Spectrum » mode then, press on F4 key [*Open*].

- If necessary, it is possible to select a different memory location for the spectrum with F1 key [*Location*] (USB medium at the USB-A connection).
- Select the required spectrum. The original view of the curve is displayed.

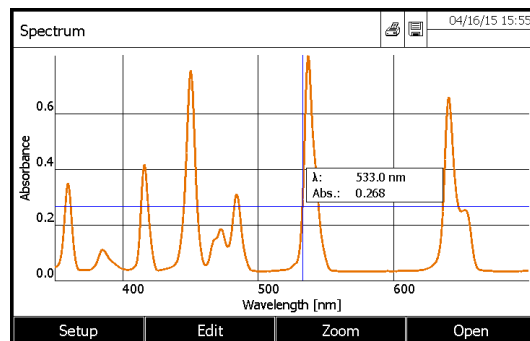1.42.2 CURSOR


The cursor consists of a horizontal and vertical line that crosses each other at a point of the curve. A box names the x and y values of the point of the curve.

Move the cursor along the x axis (wavelength) with «*◀*» and «*▶*» keys. It is possible to trace and evaluate the curve point by point.

1.42.3 ZOOM

- Press on F3 key [*Zoom*].

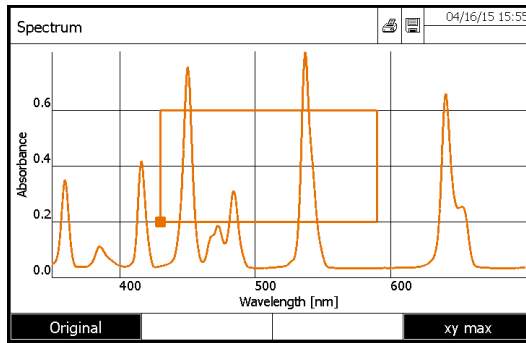UviLine Series

The zoom window appears.

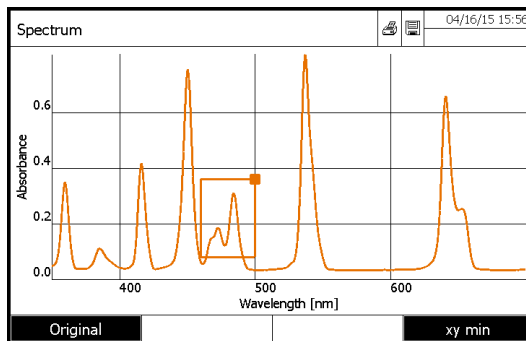
– Press F1 key [*Original*], at any time, to come back to the original view of the spectrum.

The bottom left corner of the zoom window is marked by a small black square.

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
– Use F4 [*xy max*] key to pass from the bottom left corner of the zoom window to the top right corner of the zoom window.

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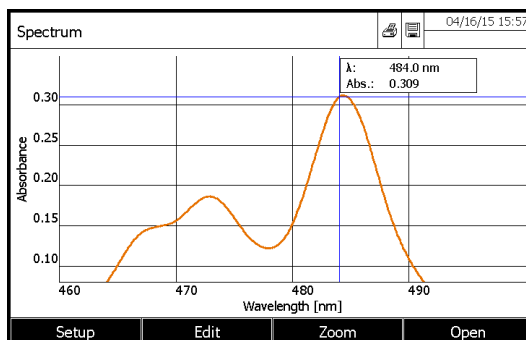– Press on <START-ENTER> key. The zoom window is scaled up on the entire diagram area.

– Press on <ESC> key to come back to the spectrum original view.

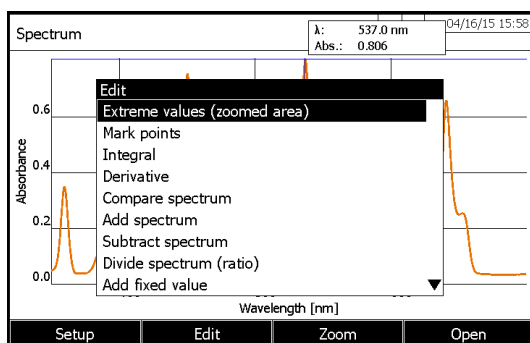
1.42.4 MANUAL CHANGE OF ABSORBANCE SCALE

Press on F1 key [Setup], select « Scaling » and then « Manual » to determine manually the absorbance scale (Y axis).
Enter manually the minimum and the maximum of absorbance.

1.42.5 CALCULATIONS

1.42.5.1 Picks and valleys detection

Press on F2 key [Edit].

Select « Extreme values » with «▲» and «▼» keys. Confirm with <START ENTER> key.
There is to way for the peak to valley detection:

1 - Automatic

Spectrum				04/16/15 18:44
Maximum values		Minimum values		
359.8 nm	0.350	351.0 nm	0.050	
385.5 nm	0.112	368.8 nm	0.038	
416.2 nm	0.417	404.6 nm	0.040	
450.4 nm	0.754	426.7 nm	0.037	
472.9 nm	0.186	461.6 nm	0.056	
484.4 nm	0.311	477.9 nm	0.122	
537.2 nm	0.809	499.7 nm	0.035	
641.1 nm	0.660	588.9 nm	0.034	
		674.1 nm	0.035	

OK Mark

2- Manual input

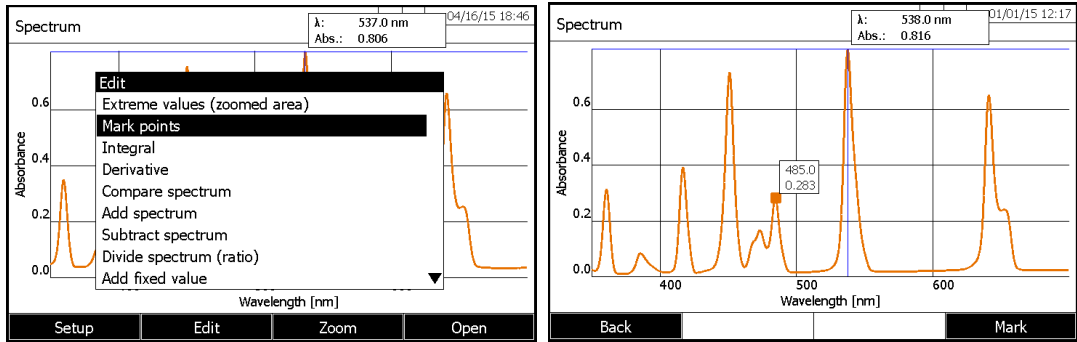
- Enter the minimum distance between minimum and maximum (More the difference between minimum and maximum will be small, more peaks and valleys will be listed)
- Press F1 key [OK] to come back to the curve.
- Press F4 key [Mark] to mark, on the curve, minimum and maximum.

1.42.5.2 Mark points


This function makes it possible to the operator to mark on the spectrum a particular point with its coordinates (Absorbance Wavelength).

- Press on F2 key [Edit].
- Select « Mark points » with «▲» and «▼» keys. Confirm with <START ENTER> key.

UviLine Series

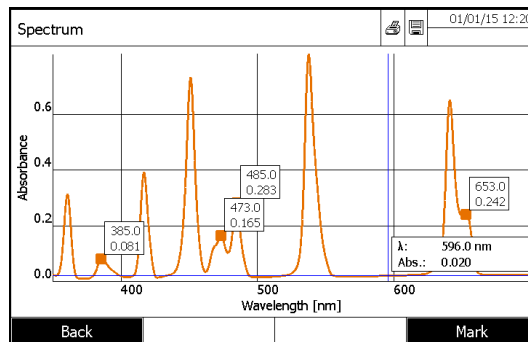
- Move the cursor along the curve with « \blacktriangleleft » and « \blacktriangleright » keys.
- Press on F4 key [Mark] to display the points coordinates.


1.42.5.3 Deleting of mark points

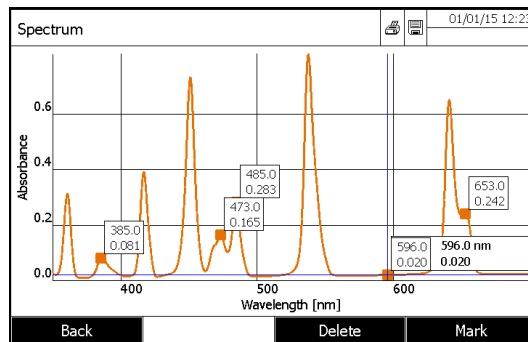Deleting of individual mark point

- Move the cursor along the curve with « \blacktriangleleft » and « \blacktriangleright » keys.
- When cursor arrives on marked point, the [Delete] key appears beside [Mark] key.
- Press F3 key [Delete] to remove the selected point.

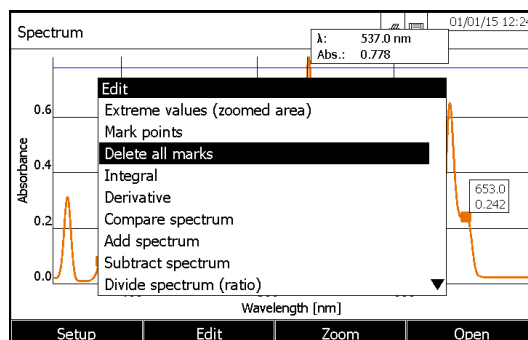Deleting of all mark points


When there are marked points, the command « *Delete all marks* » appears inside the « *Edit* » menu.

- Select « *Delete all marks* » with «▲» and «▼» keys. Confirm with <START ENTER> key. All marked points will be deleted.

1.42.5.4 Area calculation

Calculate the surface under the curve between two wavelengths [X1, X2] freely chosen by the operators.

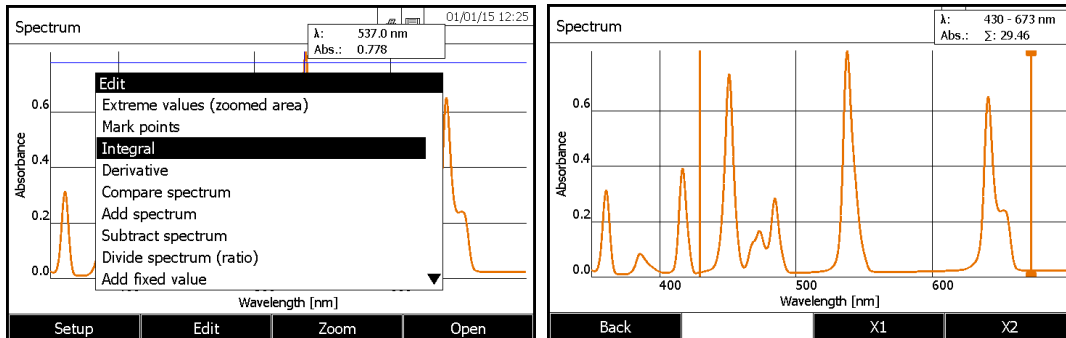- Press on F2 key [*Edit*].
- Select « *Surface calculation* » with «▲» and «▼» keys. Confirm with <START ENTER> key.


- Press on F3 key [*X1*] to select left cursor.
 - Press on F4 key [*X2*] to select right cursor.
- The surface value and the selected wavelength range are displayed at the top right corner of the screen.

1.42.5.5 Derivative calculation

Calculate the derivative of the complete spectrum. For the calculation of the second and the third derivative, it is possible to carry out the function on several times.

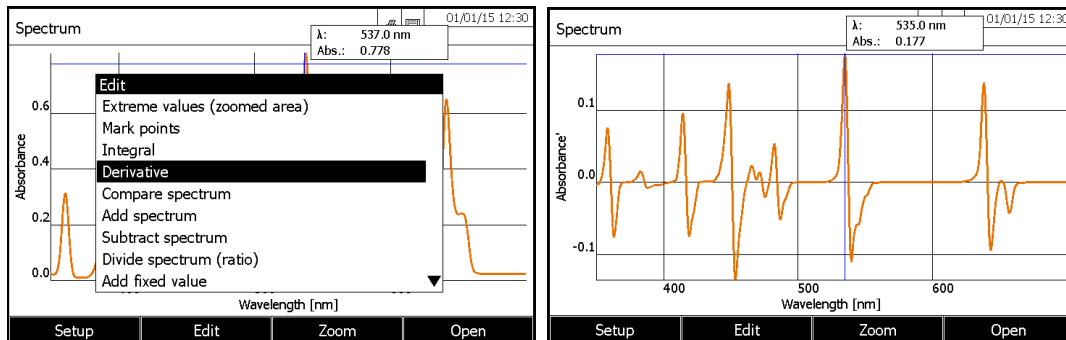
- Press on F2 key [*Edit*].
- Select « *Derivative* » with «▲» and «▼» keys. Confirm with <START ENTER> key.

Back to the original curve

- Press on F2 key [*Edit*].
- Select « *Original values* » with «▲» and «▼» keys. Confirm with <START ENTER> key.

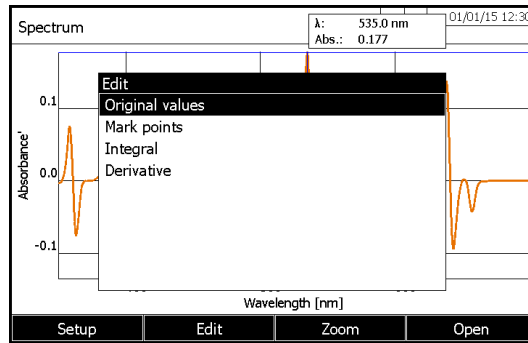UviLine Series

NOTE: The functions as « Zoom », « Mark points », « Surface calculation » and « Derivative » are also available for a derivative curve.

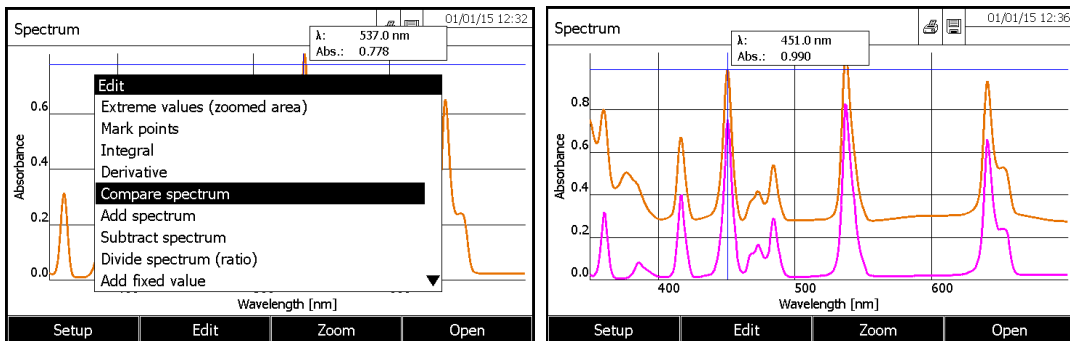
A derivative curve can be also recorded in memory with <STORE> key.

1.42.5.6 Spectra comparison

Load a second spectrum into the same diagram for direct comparison.

NOTE: The comparisons of spectra can be done only on spectra which do not result from calculations (derivative, multiplication...)

- Press F4 key [Open] and select the first spectrum
- Press F2 key [Edit].
- Select « Compare spectrum » with «▲» and «▼» keys. Confirm with <START ENTER> key.
- Select the second spectrum and confirm with <START ENTER> key.


To come back to a normal screen (first selected spectrum):

- Press on F2 key [Edit].
- Select « Original values » with «▲» and «▼» keys. Confirm with <START ENTER> key.

1.42.5.7 Addition of two spectra

- Press F4 key [Open] and select the first spectrum or use the acquired spectrum
- Press F2 key [Edit].
- Select « Add spectrum » with «▲» and «▼» keys. Confirm with <START ENTER> key.
- Select the second spectrum and confirm with <START ENTER> key.

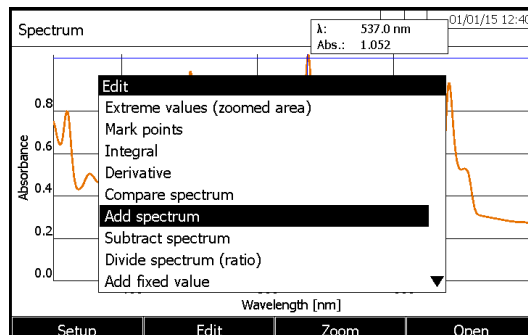It displays the summary of the two spectra.


To come back to a normal screen (first selected spectrum):

- Press on F2 key [Edit].
- Select « Original values » with «▲» and «▼» keys. Confirm with <START ENTER> key.

1.42.5.8 Subtraction of one spectrum from another one

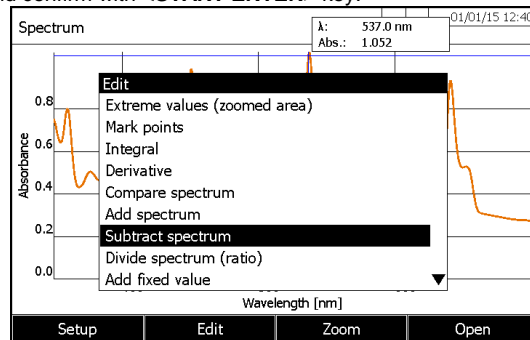- Press F4 key [Open] and select the first spectrum or use the acquired spectrum
- Press F2 key [Edit].
- Select « Subtract spectrum » with «▲» and «▼» keys. Confirm with <START ENTER> key.
- Select the second spectrum and confirm with <START ENTER> key.


To come back to a normal screen (first selected spectrum):

- Press on F2 key [Edit].
- Select « Original values » with «▲» and «▼» keys. Confirm with <START ENTER> key.

1.42.5.9 Division of a spectrum

Divide the absorbance or % transmission values of the current spectrum by the values of a stored spectrum

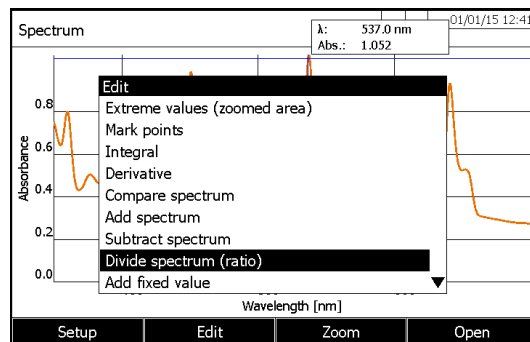- Press F4 key [Open] and select the first spectrum or use the acquired spectrum
- Press F2 key [Edit].
- Select « Divide spectrum (ratio) » with «▲» and «▼» keys. Confirm with <START ENTER> key.
- Select the second spectrum and confirm with <START ENTER> key.


To come back to a normal screen (first selected spectrum):

- Press on F2 key [Edit].
- Select « Original values » with «▲» and «▼» keys. Confirm with <START ENTER> key.

Note: This function always applies to the common wavelength range of both spectra only.

1.42.5.10 Addition of a constant

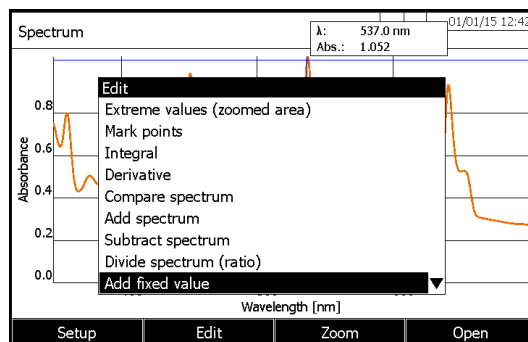- Press F2 key [Edit].
- Select « Add fixed value » with «▲» and «▼» keys. Confirm with <START ENTER> key.
- Enter the positive or negative fixed value and confirm with <START ENTER> key.


To come back to a normal screen:

- Press on F2 key [Edit].
- Select « Original values » with «▲» and «▼» keys. Confirm with <START ENTER> key.

1.42.5.11 Multiplication by a constant

Multiplies the absorbance or % transmission values of the current spectrum by a constant value.

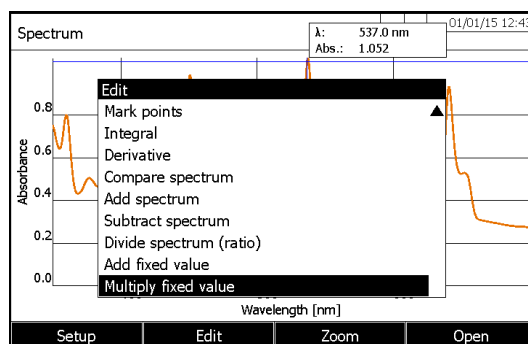- Press F4 key [Open] and select the first spectrum or use the acquired spectrum
- Press F2 key [Edit].
- Select « Multiply fixed value » with «▲» and «▼» keys. Confirm with <START ENTER> key.
- Enter the constant value and confirm with <START ENTER> key.

To come back to a normal screen:

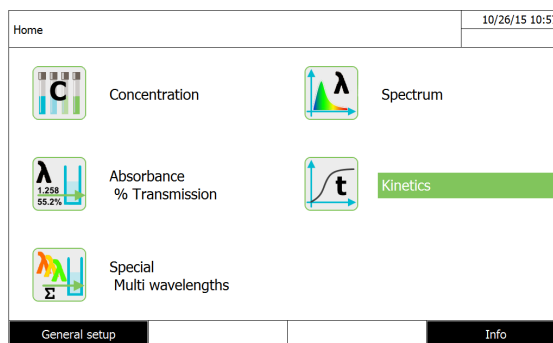
- Press on F2 key [Edit].
- Select « Original values » with «▲» and «▼» keys. Confirm with <START ENTER> key.

12 KINETICS MODE

1.43 GENERAL POINTS

The Kinetics function enables the temporal tracing of the absorbance or transmission of a sample at a certain wavelength. For recording, the spectrophotometer carries out single measurements at regular intervals (measuring interval) and stores the measured values as a time function.

All measuring parameters are entered by the operator in order to create methods which can be stored, edited and deleted.

1.44 METHOD PROGRAMMING

- Press on F1 key [New method].

Edit profile (1 of 2)		01/01/15 12:47	
Number	4000		
Name	NADH		
Mode	Absorbance		
Wavelength	525 nm		
Duration	00:10:00		
Interval	00:01:00		
Delay	00:00:00		
Colour of graph	Blue		
Scaling	Auto		
		Profile list	Delete
		Next	

Edit profile (2 of 2)		01/01/15 12:49	
Measurements/interval	1/interval		
Cell changer mode	Single cell		
Catalytic activity	Yes (only Abs)		
Factor	1.000		
Unit	µkat		
Resolution	0.01		
		Back	Complete

* **Number**
4001...4020

* **Name**
Name of the method (18 characters).

* **Mode**
Absorbance or Transmission

* **Wavelength**
320 – 1100 nm for l'UviLine 8100C / 9100C / 9300
190 – 1100 nm for l'UviLine 9400C / 9600

* **Duration (Delay + Measure)**
Total duration of the kinetics including the waiting time (Delay).
Format hh:mm:ss (hours:minutes:seconds)

* **Interval**
It is the time interval between two successive single measurements.
Format hh:mm:ss (hours:minutes:seconds)

* **Delay**
Time between the start of the recording and the start of the first single measurement. This time allows avoiding the no linear part of a reaction. Format hh:mm:ss (hours:minutes:seconds)

*** Scaling**

Auto or Manual

*Scaling: Auto***

The spectrophotometer adjusts the scaling of the axes (minimum and maximum value of the axis) to the measured values while measuring. The entire curve is always visible.

Scaling: Manual

Abs min.

Abs max.

The scaling of the axes (minimum and maximum value of the axis) is permanently set manually.

*** Measurements/interval**

1/interval or Max./interval

Here, it is possible to define how many measurements will be carried out per interval. The setting affects the calculation of the slope of the individual intervals.

The displayed slope for an interval is determined as follows, depending on the profile:

Measurements/interval	Slopes
1/interval	Difference of two measured values (2 points calculation)
Max./interval	Slope of the straight line determined in an interval by linear regression

*** Katalytic activity**

YES or NO

Katalytic activity	Result
NO	Total slope report to the minute
YES	Total slope report to the minute multiplied by a factor

*** Factor**

Factor multiplying the average slope report to the minute.

*** Units**

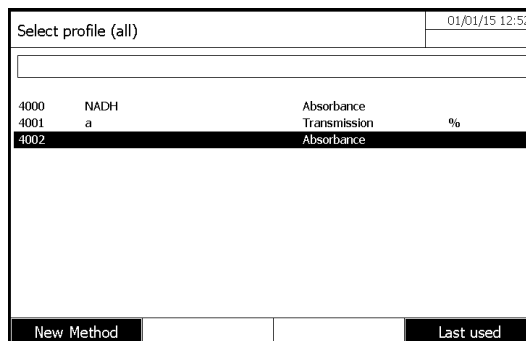
i.e.: μkat (10 characters maximum).

*** Resolution**

From 0 to 3

Number of digits, after comma, for the result.

1.45 METHOD SELECTION

- Select a method in the "Kinetics" mode then press on **<START-ENTER>** key to confirm the method choice.

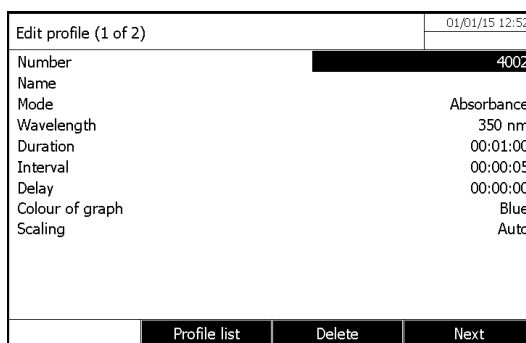
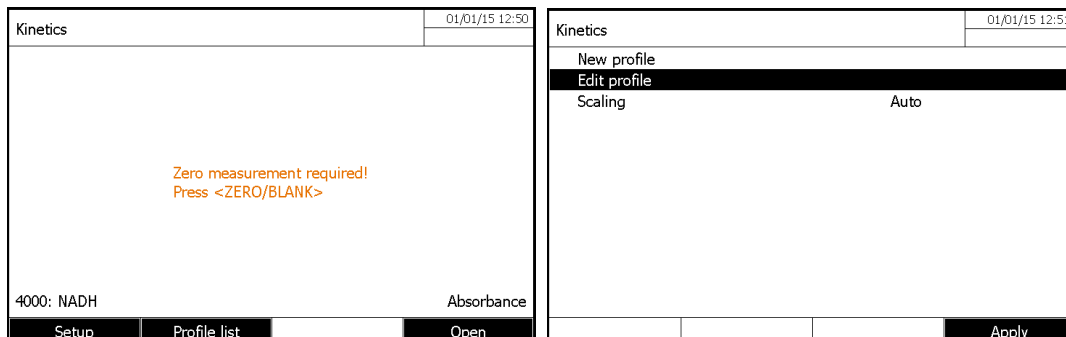
1.46 METHOD SAVING

Methods are automatically saved in memory.

1.47 METHOD DELETING

- Select the method to be deleted with «▲» and «▼» keys. Confirm with <START ENTER> key. Then press F1 key [Setup].

- Select "Edit method" with «▲» and «▼» keys. Confirm with <START ENTER> key.

- Press F3 key [Delete].

1.48 LIMITATION OF METHODS LIST

If the list is very long, it is possible to narrow down the methods list and thus make the search easier as follows:

- Press on F4 key [Last used]; it is possible to restrict the method list to the ten methods last used.

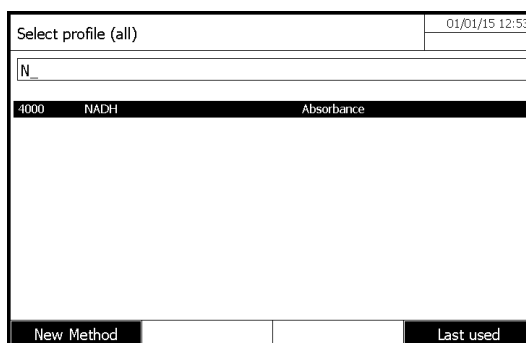
- With the search function, it is possible to search certain character strings in the list. The search takes place as a full-text search of the entire list contents.

Thus the operator can search for a method number or name.

Search for a character string:


Enter the character string to be searched for in the search window with <A...9>.

The list appearing below shows all methods containing the character string. This methods list is updated with each character that is entered.

1.49 METHOD EXECUTION

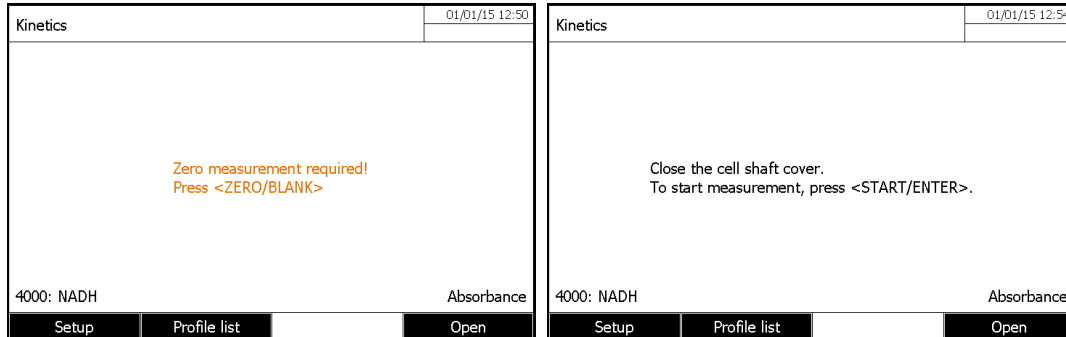In the "Kinetics mode" select a method with «▲» and «▼» keys and- Press on <START-ENTER> key to confirm the method choice. Then -press <ZERO BLANK> key


- Insert sample.
- Press <START ENTER> key to start the measurement. The spectrophotometer starts recording automatically.
- Wait the end of measurement.

Stopping the recording:

- Press on F4 key [*Stop*] to terminate the recording prematurely. The curve recorded up to this point can be stored and edited (see paragraph 12.8.3).
- Press on <ESC> key to completely cancel measurement. The curve recorded up to this point is discarded.

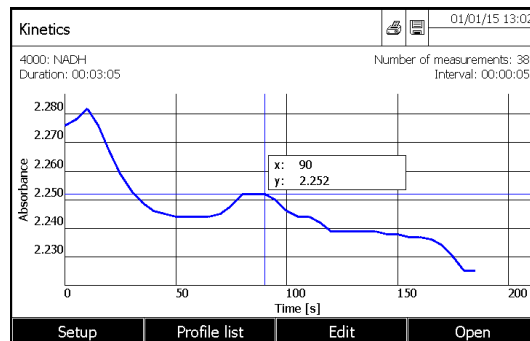After the specified *Duration* has expired, the cursor appears.

- Press on <START ENTER> key to run a new measurement.

1.50 METHOD EDITING

- Enter in « Kinetics » mode.
- Select a method with «▲» and «▼» keys, confirm with <START ENTER> key.
- Press F1 key [Setup].
- Select « Edit method » with «▲» and «▼» keys, confirm with <START ENTER> key.

1.51 KINETICS TREATMENT

1.51.1 KINETIC CURVE PRINTING

- Press <PRINT> key to output the kinetic curve to a connected printer as a graphic.

1.51.2 CURSOR

The cursor consists of a horizontal and vertical line that crosses each other on a point of the curve. A box names the x and y values of the point of the curve.

Move the cursor along the x axis (time axis) with «◀» and «▶» keys. By this way, it is possible to trace and evaluate the curve point by point.

These points are corresponding to the measurement interval.

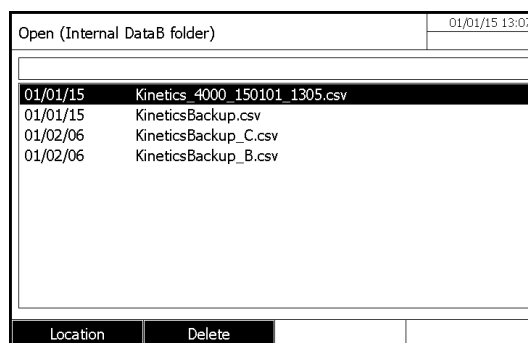
1.51.3 KINETIC CURVE SAVING

Press <STORE> key to save the kinetic curve.

If the operator forgot to save the curve, this one is nevertheless kept in memory under the name of "date kineticsBackup.csv".

1.51.4 DOWNLOADING OF STORED KINETIC CURVE


- Press F4 key [Open].

The list with the stored kinetic records (*Internal DataB folder*) is displayed.

- Press F1 key [Location] to select another memory location of the kinetic record (*USB memory*)
- Select the kinetic curve and validate.

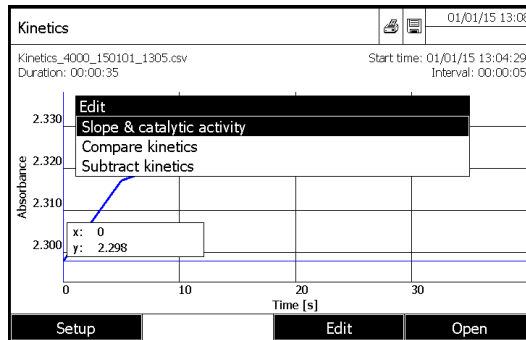1.51.5 « EDIT » MENU

1.51.5.1 Slope

Select « Slope & catalytic activity ».

Interval	Slope [Δ/min]	Time [s]
1	0.228	5 s
2	0.060	10 s
3	0.024	15 s
4	0.036	20 s
5	0.036	25 s
6	0.012	30 s
7	0.024	35 s
Total	0.049	35 s

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It is the number of measurement interval.

« Slope [Δ/min] » column

It is the slope report to the minute per interval.

« Time [s] » column

It is the time at the end of each interval.

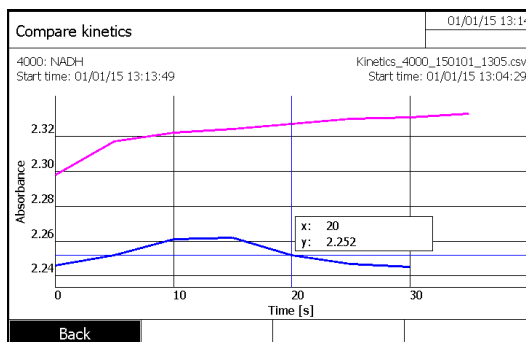
RESULT:

It is the slope, report to the minute, of the straight regression line passing through various points of measurement, multiplied by the factor and expressed with the resolution and unit selected by the operator.

1.51.5.2 Comparison of kinetic curves

Load a second kinetics curve into the same diagram for direct comparison.

- Press on F3 key [Edit].
- Select « Compare kinetics ».
- Select another kinetics curve. Confirm with <START ENTER> key

To come back to a normal screen (first selected kinetics curve):
- Press on F1 key [Back].

1.51.5.3 Subtraction of a kinetic curve from another one

Subtract a stored kinetics curve from the current kinetics curve.

- Press on F3 key [Edit]
 - Select « Subtract kinetics ».
 - Select a kinetics curve. Confirm with <START ENTER> key.
- The curve resulting of this operation can't be saved in memory.
To come back to a normal screen (first selected kinetics curve):
- Press on F1 key [Back].

NOTE: To subtract one kinetics curve from another one, the intervals have to be identical.

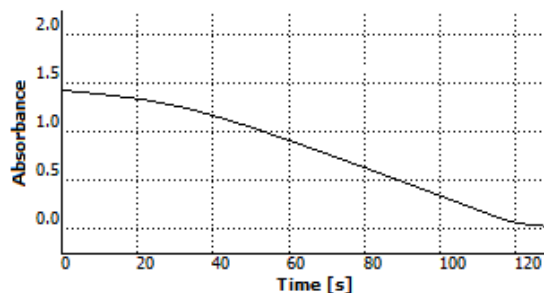
1.51.6 MANUAL CHANGE OF ABSORBANCE SCALE

Press on F1 key [Setup], select « Scaling » and then « Manual » to determine manually the absorbances scales (Y axis).
Enter the absorbance minimum and maximum.

1.52 CURVE PRINTING


- Press <PRINT> key to print the curve on a connected printer.

UviLine 9300 7449001 1.14
Administrator
04/10/2015 14:05:46
4003

13 MEMORY

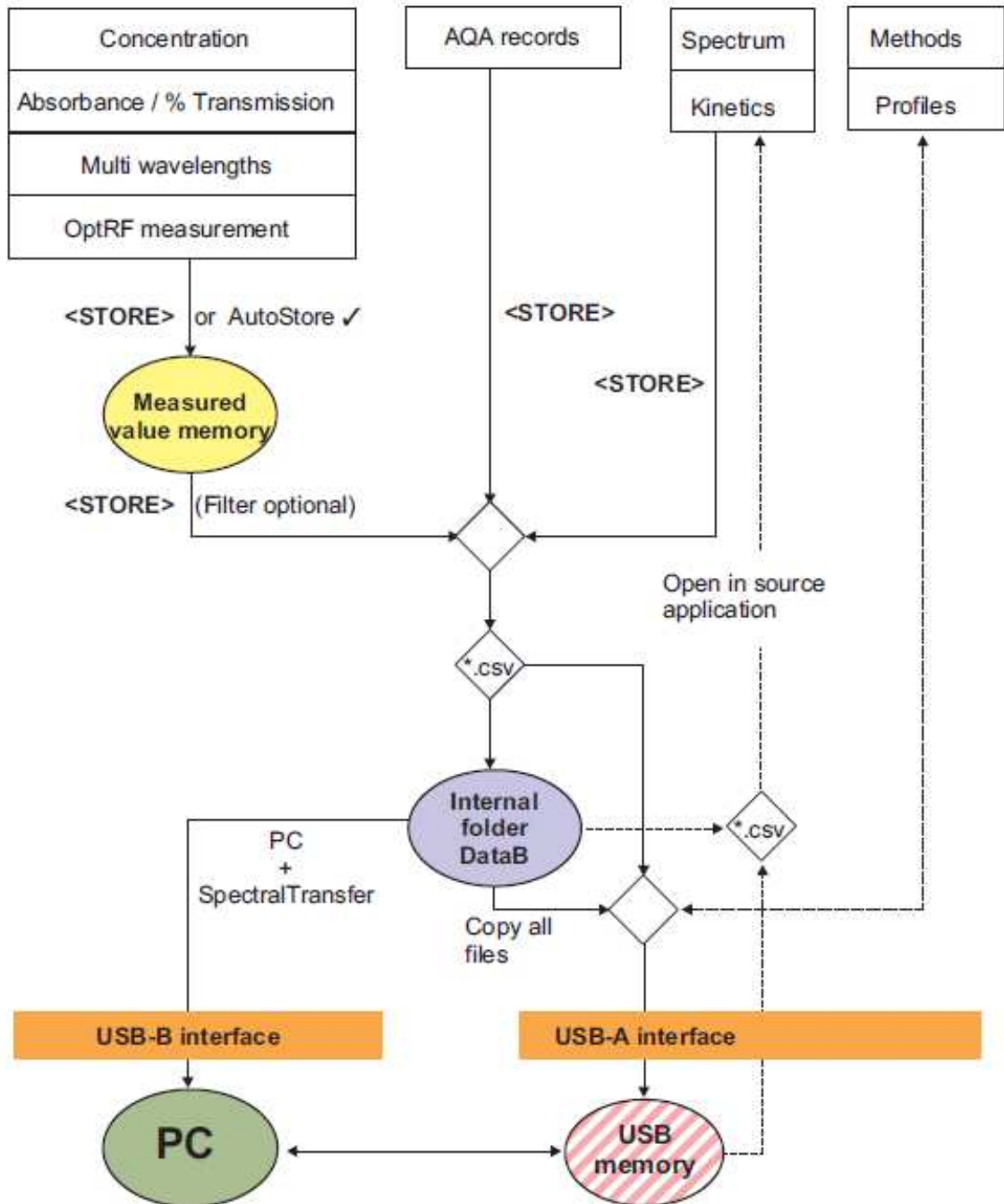1.53 OVERVIEW

Measurement data	Saving and exporting
Concentration Absorption / % transmission Multi-wavelengths	<p>Measurement datasets of these modes are saved with <STORE> or <i>AutoStore</i> in the measured value memory of the spectrophotometer first (5000 locations).</p> <p>The measured value memory is available via the <i>Measurement data memory</i> menu.</p> <p>Here it is possible to view or filter stored measurement datasets or export them into a PC readable file (*.csv) (<STORE>).</p> <p>Csv files of these modes cannot be imported back into the photometer.</p>
Spectrum Kinetics	<p>Measurement datasets of these modes are saved and exported as a PC readable file (*.csv) with <STORE>.</p> <p>Csv-files of these modes can be re-imported and viewed on the photometer.</p>

As the location for the PC readable files (*.csv) to be stored, it is possible to select for each export either the spectrophotometer (*Internal DataB folder*) or an external memory (*USB memory*).

Later it is possible to transmit the files stored in the photometer (*Internal DataB folder*) to a connected PC or an external memory (*USB memory*).

It is possible to transmit to a USB memory either individual files with measuring data or all files stored in the *Internal DataB folder*.

1.54 INSTRUCTIONS ON USING USB MEMORY DEVICES

The safety of data stored on USB memory devices depends on the quality of the memory device and the data transmission.

Data is stored partly or not at all if for example:

- The power supply of the external memory device is interrupted during the write process, or
- The external memory device is prematurely disconnected from the photometer during the data backup.

To prevent a data loss we recommend the following:

- Save all data internally in the photometer first.
- After performing a backup leave the USB memory device connected to the photometer for some time.
- Check whether the stored data is complete, e.g. on a PC.
- Use the USB memory device for data transport but not for permanent data storage.

1.55 MEASUREMENT DATASETS

1.55.1 ELEMENT OF A MEASUREMENT DATASET

A complete measurement dataset consists of:

- Consecutive number (is automatically assigned by the spectrophotometer).
- Date/time.
- Identification (e.g. ID or "AutoStore").
- User name.
- Measured parameter, e.g. method number, dilution, wavelength (depending on the measuring mode).
- Measured value with unit and if necessary citation form.

1.55.2 OPÉRATIONS WITH MEASUREMENT DATASETS

Measurement datasets can be

- Stored
- Displayed and printed
- Filtered, i.e. selected or hidden based on certain criteria
- Deleted
- Copied

UviLine Series

Remark:

If the memory is full, you can erase measurement datasets, or overwrite the oldest dataset with the next storing procedure. A security prompt appears before a dataset is overwritten. To backup the measurement data, you can transmit the measurement datasets from the measurement data memory to the internal DataB folder or a USB memory device connected to the USB-A connection and archive them further from there

1.55.3 RESULT RECORDING

1.55.3.1 Manual recording

After each measurement, it is possible to store the measurement data manually with the <STORE> key. It is stored in the measurement data memory. The memory symbol  in the header indicates that the measurement data displayed on the screen is ready to be stored. With the measuring modes, *Concentration*, *Absorbance / % Transmission*, and *Multi wavelengths* there is the additional option to automatically store all new measured values at the time of the measurement

When storing manually, an input field for the identification (ID) appears after pressing the <STORE> key. Here it is possible to enter an individual combination of alphanumeric characters for later easier identification of the measurement datasets. 30 digits are available for this.

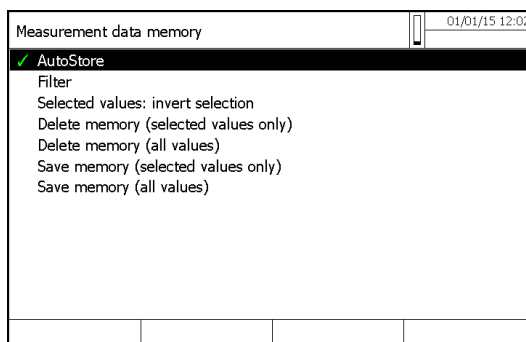
WARNING:

In « Absorption / % transmission » mode, the « AutoStore » function is not operational for data acquired during this measurement.

If the operator wants to store data, he has to use <STORE> key. So, all results acquired during last measurement will be stored under the names given by the operator.

1.55.3.2 Automatic recording with « AutoStore » function

- Press on F1 key [Setup].
- Select « AutoStore » with «▲» and «▼» keys and confirm with <START ENTER> key.

The symbol ✓ at the end of « AutoStore » indicates that the function is active. If this function is not active, it means that data will be saved manually only with <STORE> key.

Remark

For the measuring modes, *Concentration*, *Absorbance / % Transmission*, and *Multi wavelengths*, it is possible to save every measured value automatically with (AutoStore) function.

All automatically saved measurement datasets are given the label "AutoStore". If the same measurement dataset is manually saved using (<STORE>) key, the "AutoStore" label is overwritten by ID code given by the operator.

The AutoStore setting is valid for all three measuring modes, *Concentration*, *Absorbance / % Transmission*, and *Multi wavelengths*.

1.55.4 RESULT DISPLAYING

From main menu

<HOME>
- [General setup]
- Measured value memory

From concentration menu

<HOME>
- Concentration
- Choice of a memorized method
- [Setup]
- Measured value memory

From Absorption/Transmission menu

<HOME>
- Absorption/Transmission
- [Setup]
- Measured value memory

From Multi-wavelengths menu

<HOME>
- Multi-wavelengths
- Choice of a memorized method
- [Setup]
- Measured value memory

Each one of these possibilities displays the contents of the measurement data memory in the form of list in the following way:

Concentration	01/01/15 13:15
Dilution	
Sample blank value	
✓ Display absorbance	
New method	
Edit method	
Delete method	
Measurement data memory	

- Select « Measurement data memory » with «▲» and «▼» keys and confirm with <START ENTER> key.

UviLine Series

Measurement data memory

Measurement date & time	Result of the measurement	Citation form
01/01/15 12:09	20.93 mg/l NO3	Test2
01/01/15 12:08	> 100.00 mg/l NO3	AutoStore
01/01/15 12:08	> 100.00 mg/l NO3	Test1
01/01/15 12:08	> 100.00 mg/l NO3	AutoStore
01/01/15 14:47	-0.000 Absorbance	Sample 5
01/01/15 14:47	0.001 Absorbance	Sample 4
01/01/15 14:47	-0.006 Absorbance	Sample 3
01/01/15 14:47	0.001 Absorbance	Sample 2
01/01/15 14:47	0.001 Absorbance	Sample 1
01/01/15 12:03	0.001 Absorbance	AutoStore
01/01/15 12:18	0.000 Absorbance	AutoStore
01/01/15 12:18	-0.000 Absorbance	AutoStore
01/01/15 12:17	0.000 Absorbance	AutoStore
01/01/15 12:17	0.009 Absorbance	AutoStore
01/01/15 12:17	-0.000 Absorbance	AutoStore

Memory space usage: 369/5000

Buttons: Setup, Single value, Delete, Save

If there are more datasets available than can be displayed, the arrows «▲» and «▼» displayed additionally.

Filter ✓ indicates that filter settings are active. In this case, only those datasets are displayed that correspond to the selected filter criteria.

Options:

Measurement datasets can be:

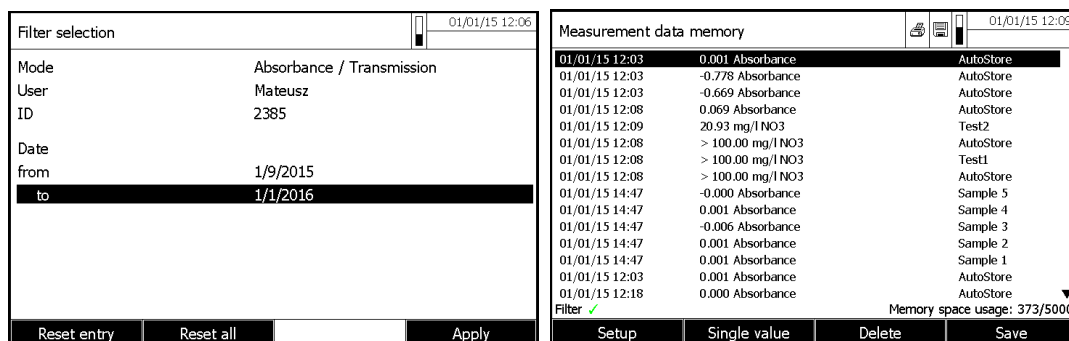
- displayed in short form as a list or in details as individual values ([List] <-> [Single value])
- filtered (see section 4.12.7 and section 4.12.8)
- deleted (see section 4.12.9).
- with <STORE>, you can store the entire displayed list as a *.csv file in the internal DataB folder or on a USB memory device connected to the USBA connection. The filter settings apply to the storing process. You can freely select the file name. Thus you can, e. g. store in a separate file and systematically archive measurement data of a certain period.
- with <PRINT>, the entire displayed list can be printed. The filter settings apply to the print process.

1.55.5 RESULT FILTERING

1.55.5.1 Filtering of a results group

It is possible to select a result or a group of results with the "Filter" function.

The functions display, delete and download stored measurement datasets refer to all stored measurement datasets that correspond to the specified filter criteria.

Filter criteria:

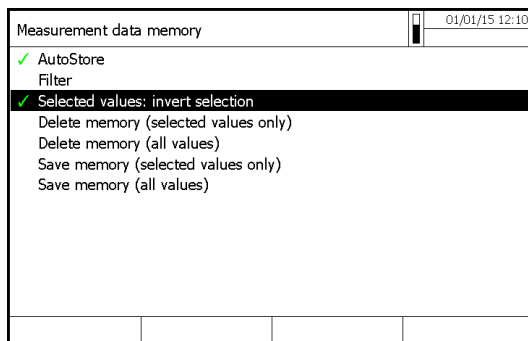
The following filter criteria can be set:

- *Mode (Absorption / % transmission and Multi-wavelengths)*
- *User*
- *ID (identification)*
- *Date (date from ... to ...)*
- *Method (for the measured parameters, Concentration and Multi-wavelength)*

Remark: Alternatively, it is possible to hide measurement datasets that meet the specified filter criteria with the "Selected values: invert selection" function).

1.55.5.2 Filters inversion

With the "Selected values: invert selection" function, it is possible to hide all measurement datasets that correspond to the specified criteria of the filter.


Remark: It is possible to use this function to select and delete measurement datasets no longer used.

1.55.5.3 Erase stored measurement dataset

It will delete only results corresponding to the normal or invert selection criteria.

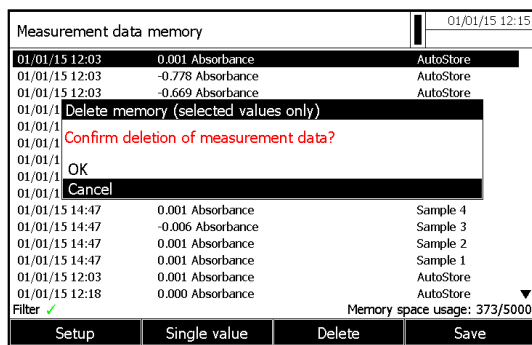
- Select « Delete memory (selected values only or all values) » with «▲» and «▼» keys and confirm with <START ENTER> key.
- Select « OK » or « Cancel »

1.55.5.4 Result deleting

It is possible to delete only one result.

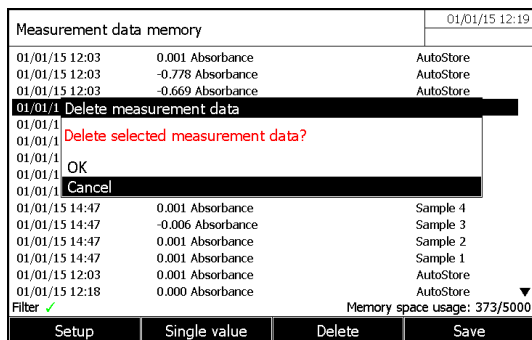- Select the result to be deleted with «▲» and «▼» keys in the data memory.
- Press F3 key [Delete].
- Select « OK » or « Cancel »

1.55.5.5 Saving Kinetic recordings, spectrUM

After the following measurements, the Save dialog opens and prompts you to save the data in a *.csv file:

- Kinetics
- Spectrum

If the data are not saved in *.csv format, they are lost when the measuring mode is terminated.

Remark:

During a kinetic recording, the current measurement is always saved in the file, "KineticsBackup.csv" for safety reasons.

1.55.5.6 Saving data as a pdf file

All data that can be printed (printer symbol on the display) can also be saved as a pdf file. The pdf file contains the data that are also output to a USB printer. Kinetic recordings and spectra are stored in the pdf file as a graphic.

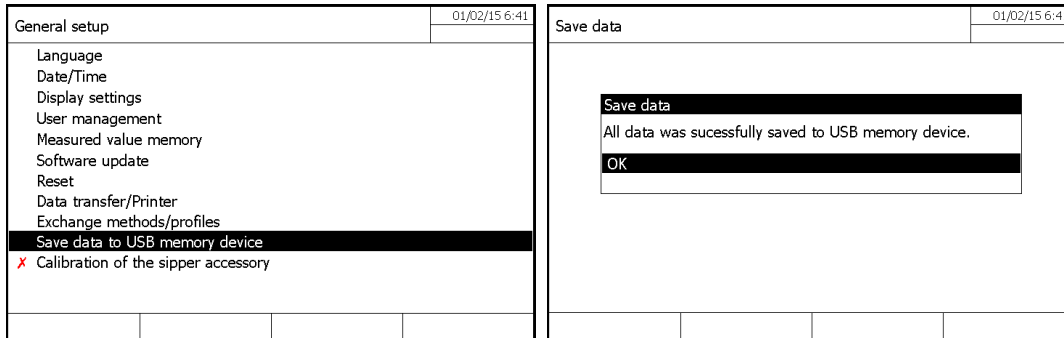
Saving as a pdf file and printing is done with the <PRINT> key. Prerequisite is that pdf printing is set as the printer in the menu <HOME>/General setup/ Data transfer/Printer/Function of PRINT key. Subsequently, enter a file name and select the storage location (internal folder DataB or USB memory device).

1.55.6 SAVING / EXPORTING FILES

If you want to back up or process measurement data files outside the photometer, you can copy them to external media.


1.55.6.1 Copying all measurement data files to a USB memory device

You can very simply transfer all measurement data files from the photometer (Internal DataB folder) to a connected USB memory device.

All measurement data files from the photometer (Internal DataB folder) have been transferred to the USB memory device. The complete folder structure from the photometer is created on the USB memory device. The individual measurement data files are stored in subfolders sorted by measurement data types.

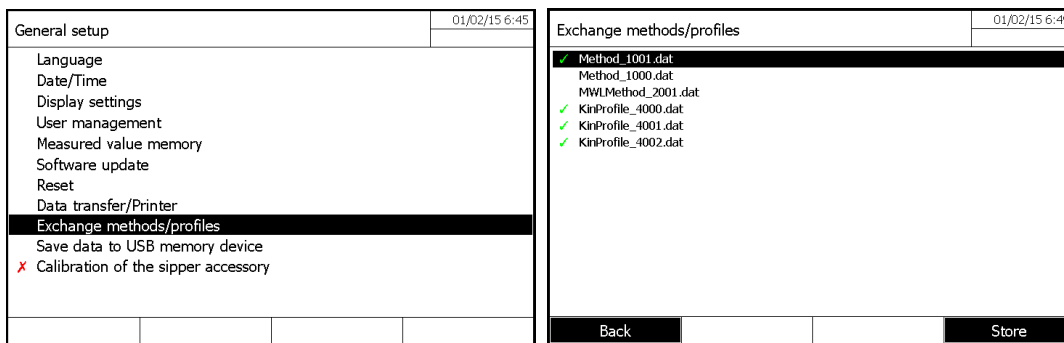1.55.6.2 Copying user-defined methods / profiles to a USB memory device

A list is displayed that includes all user-defined methods and profiles available on the photometer. All methods and profiles are checked off with a checkmark. All methods and profiles checked off are saved.

If necessary, select individual methods/profiles with and remove the checkmark with **<STAR/ENTER>**. These methods/profiles will not be saved. Start the save process with **<STORE>**.

A message appears when the data have been saved. Confirm the message with **<START/ENTER>**.

The save process is completed. The data are stored in the Exchange_Method_Profile folder on the USB memory device. The individual files with the methods/ profiles are in subfolders. Already existing files with identical names are overwritten without confirmation prompt

1.55.6.3 Copy of individual file on USB memory

1.1.1.1.1 Spectrum and kinetic curve after acquisition

After measurements of spectrum or kinetics, the user can save his curve spectral or kinetic in a "*.csv" file:

- Press F4 key [*Open*] to enter in the safeguard sub-menu.
- Press F1 key [*Location*].
- Select « USB memory ».
- Give a name.
- Validate with **<START/ENTER>** key.

If the safeguard, in "*.csv" format were not carried out, the measurement data are lost when the user leaves the measurement mode.

UviLine Series

Remark

With kinetic recordings, the current measurement is always saved in the "KineticsBackup.csv" file for safety reasons.

1.1.1.1.2 Spectrum and kinetic curve already memorized inside the « Internal file DataB »

- Press F4 key [Open].
- Select a spectrum or kinetics curve.
- Confirm with <START/ENTER> key.
- Press <STORE> key to open save dialog box.
- Press F1 key [Location].
- Select « USB memory ».
- Change the name if necessary.
- Validate with <START ENTER> key.

1.1.1.1.3 Absorption/Transmission, Concentration, Multi-wavelengths mode

The measurement data of Absorbance/ Transmission, Concentration and multi-wavelengths are stored in the measurement data memory first, either automatically or manually (with the <STORE> key). Data stored in the measurement data memory can be filtered with filter criteria and then exported in the PC readable in "*.csv" format.

- Press F1 key [Setup] inside one of the measuring mode or [General setup] from the main menu
- If necessary, set any filter criteria with [Setup].
- Press <STORE> key to open a save dialog box.

The spectrophotometer automatically suggests the location, *Internal DataB folder*, and a file name.

- If necessary, use F1 key [Location] to change the location (USB memory).
- If necessary, change the suggested file name.
- Press <START ENTER> to save the selected results.

1.55.7 PRINTING THE DATA (USB)

Usable printers

Data can be printed with standard printers (ink-jet or laser) connected to the USB-A interface. Suitable are all printers that can interpret the PCL-3 printer control language.

The printer symbol  indicates that the display contents can be printed. To print, press <PRINT>.

Pdf file

As an alternative, you can also output the print data to a pdf file.

Remark

In den following paragraphs, "Print" means:

- output to a USB printer
- output to a pdf file.

1.55.8 SETTING FOR DATA TRANSMISSION

Decimal separator for csv-Files

For the output of CSV files, it is possible to select a comma or point as the decimal separator.

Data format

When printing measurement datasets, it is possible to select a short or long version with different information contents.

Function of PRINT Key

Her you can choose between a USB printer & PDF file

1.55.9 IMPORTING FILES

You can import to an UviLine spectrophotometer the data that were created with the same or another UviLine spectrophotometer, and were then saved to a USB memory device.

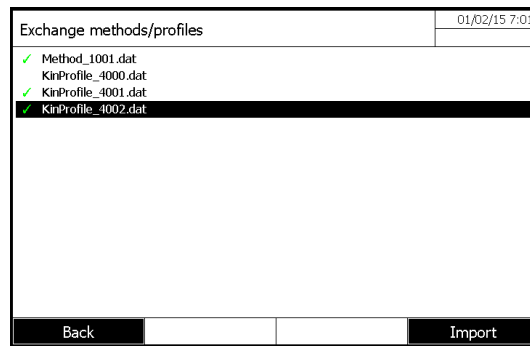
You can import the following data:

- Spectra
- Kinetic recordings
- User-defined methods
- Profiles

1.55.9.1 Importing spectra or kinetic recordings from a USB memory device

You can import to the photometer any spectrum or kinetic recording by opening an externally stored spectrum or kinetic recording with the Open function of the photometer.

1.55.9.2 Importing methods / profiles from a USB memory device

A list is displayed including all user-defined methods and profiles stored in the corresponding subfolders of the Exchange directory on the USB memory device. All methods and profiles are checked off with a checkmark. All methods and profiles checked off are imported.

Start the import with [Import]. A confirmation prompt appears before any data on the photometer are overwritten. A message appears when the data have been imported.

Confirm the message with <START-ENTER>. The import is completed. The imported methods / profiles are available on the photometer.

Remark

When importing methods make sure that your photometer supports the wavelengths of the imported methods.

14 USERS MANAGEMENT

There are a user name and a password by fault:

- User name: **Administrator**
- Password: **admin**

Respect the spaces and the letters size (Capital or small letter).

The functions of the user management are only available for users of the user group, *Administrator*.
An administrator can:

- Activate or deactivate the user management for the spectrophotometer.
- Create, change or delete individual user accounts.

The spectrophotometers UviLine Series allow the management of up to 100 users. Each user is a member of a user group with defined user rights.

1.56 USER LEVEL AND USER RIGHTS

There are three hierarchical user groups:

- *Administrator* (top level)
- *User* (user account registered by the administrator)
- *Guest* (user without user account)

Administrators and users log in to the photometer with their user name and password.

Guests can optionally enter a name for their login. Thus, documented measured values can later be assigned to the user.

Action	Administrator	User	Guest
Select methods	✓	✓	✓
Execution of methods	✓	✓	✓
Record of results	✓	✓	✓
Deactivate automatic records of the results	✓	⊘	⊘
Activate or deactivate « users management » function	✓	⊘	⊘
Edit and modify a method stored in memory	✓	⊘	⊘
Delete a method stored in memory	✓	⊘	⊘
Date and time setting	✓	⊘	⊘
Delete of results stored in memory	✓	⊘	⊘
Users management	✓	⊘	⊘
Reset of spectrophotometer setting	✓	⊘	⊘
Update of the software	✓	⊘	⊘

Remark

You can also switch off the user management and reactivate it as necessary. To do so, you need administrator rights. If the user management is switched off, the user name and password do not have to be entered. Each user has full rights.

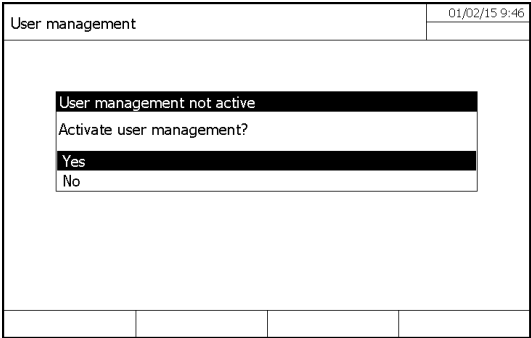
1.57 ACTIVATION/DEACTIVATION OF USER MANAGEMENT

Each user can activate the user management function. If the function is deactivated, each user has administrator rights. Only members of the user group, administrator can deactivate the user management function.

If the function is active, each user has to log in the spectrophotometer. After the login, the user has certain rights depending on the user group.


1.57.1 ACTIVATION OF USER MANAGEMENT FUNCTION

User management activation is accessible in the "general setup" menu

1.57.2 DESACTIVATION OF USER MANAGEMENT FUNCTION

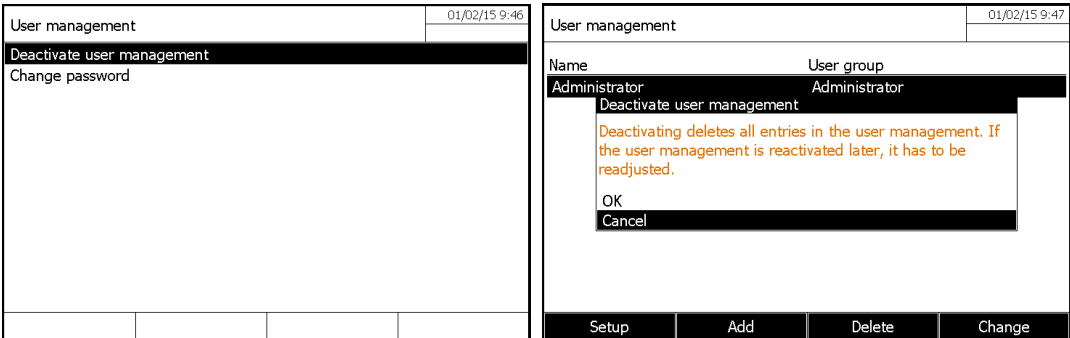Select « Deactivate user management » and confirm. Then, select « OK » and confirm. The user management function is deactivated. Each user has administrator rights.

REMARK

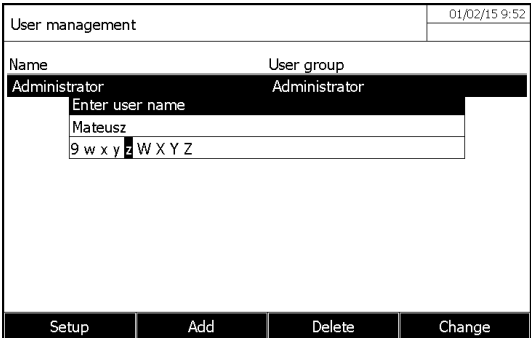
If the user management is deactivated by a user of the *Administrator* user group, all user accounts that were set up are lost. The password for the administrator is reset to "admin".

1.58 CREATION, MODIFICATION OR DELETING A USER ACCOUNT

When the user management function is active, a user with administrator rights can manage user accounts.


1.58.1 USER ACCOUNT CREATION

During the creation of a user account, the *Name*, *User group* and *Password* are defined.

1.58.2 USER ACCOUNT MODIFICATION

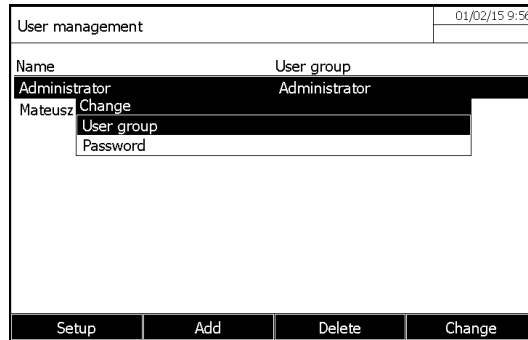At the time of the modification of a user account, it is possible to modify the membership of *User group* and the *Password*. Only the user having an Administrator account is authorized to make these changes.


The screenshot shows a terminal window titled "User management" with a timestamp of "01/02/15 9:56". The interface displays a table with two columns: "Name" and "User group". The first row shows "Administrator" in both columns. Below the table, the name "Mateusz" is visible, and a "Change" menu is open, showing options for "User group" and "Password". At the bottom of the window, there are four buttons: "Setup", "Add", "Delete", and "Change".

1.58.3 PASSWORD MODIFICATION

The administrator creates user accounts and assigns a password on each user account. As soon as a user opened a session successfully with his user account, it can modify itself the password of his user account.

1.58.4 USER ACCOUNT SUPPRESSION

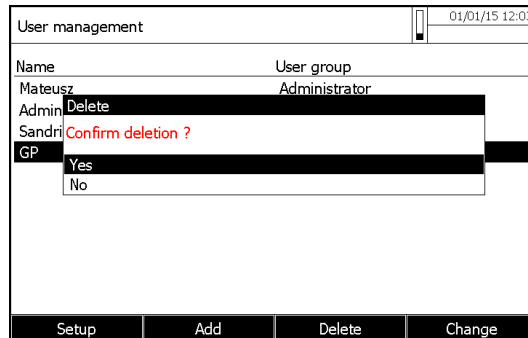1. Select a user account.
2. Press F3 key [*Delete*] to delete the user account: A security prompt appears.
3. Confirm the security prompt: The user account is deleted.

The screenshot shows a terminal window titled "User management" with a timestamp of "01/01/15 12:03". The interface displays a table with two columns: "Name" and "User group". The first row shows "Admin" in the "Name" column and "Administrator" in the "User group" column. Below the table, the name "Sandri" is visible, and a "Delete" menu is open, showing a security prompt: "Confirm deletion ?". Below the prompt, there are two options: "Yes" and "No". At the bottom of the window, there are four buttons: "Setup", "Add", "Delete", and "Change".

15 VARIOUS ADJUSTMENT

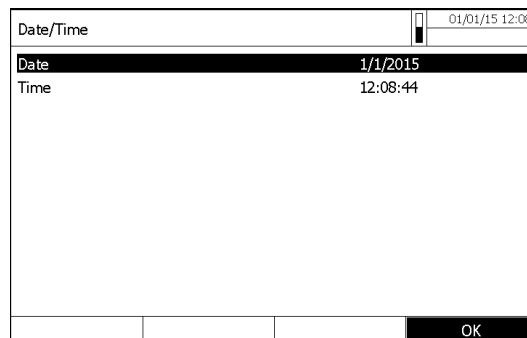
1.59 LANGUAGE CHOICE

In "general setup" menu, select "Language", then validate the selected language with <START-ENTER> key.

The language selected is noted with the symbol ✓.

1.60 DATE AND TIME ADJUSTMENT

In "general setup" menu, select "Date/Time".

Enter the current date with <0...9> keys and confirm. The input field closes. The date is accepted.

Enter the current hour with <0...9> keys and confirm. The input field closes. The time is accepted.

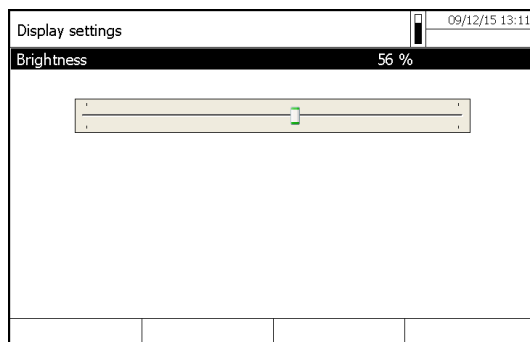
1.61 CONTRAST ADJUSTMENT

It is possible to adapt the contrast to the lighting conditions.

Select and confirm « Contrast ». An adjustment system using a cursor is displayed to adjust the screen contrast.


Adjust the screen contrast by pressing « ◀ » and « ▶ » keys and confirm with <START/ENTER> key.

Come back to the main menu with <ESC> key.

1.62 RESET

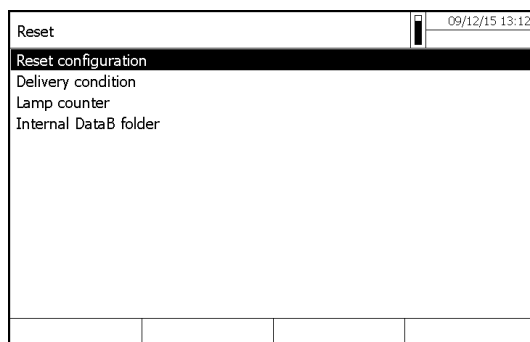It is possible to reset (initialize) the measurement settings or all settings.

Remark

The reset function is only available for users belonging to the user group « Administrator ».

There are different possibilities to reset the instruments settings.

❖ <i>Parameters reset</i>	All settings are deleted except the memory which contains measurements results, methods defined by the user and the measured blank values.
❖ <i>Delivery conditions</i>	All settings (including measurement data memory and user-defined methods) are deleted and the spectrophotometer is reset to the delivery condition.
❖ <i>Service lamp counter</i>	Allow to reset to zero the service lamp counter after the lamp changing

1.63 LAMP SERVICING COUNT DOWN

The photometer counts the operating hours of the lamp. The information on the operating hours of the lamp is given in the Info menu. The number quoted there corresponds to the number of measurements. After the lamp was exchanged you should reset the lamp counter (previous section).

1.64 SPECTROPHOTOMETER INFORMATION ([INFO])

Following information relating to the spectrophotometers are displayed by pressing F4 "info" on the main menu:

- Instrument designation.
- Instrument serial number.
- Number of the instrument software version.
- Updating date of the software.
- Hardware version
- Hardware status (for servicing).
- Number of lamp working hour.
- Results of the tests of the most important parts of the spectrophotometer.
- Memory status.

- Registered user.

Info		09/12/15 13:16
Model designation:	UviLine 9300	
Serial number:	150430007	
Software version:	2.18	
Build:	05.04.15 13:29	
Hardware version:	2-1-1/40-36/6.S77-S1+	
Hardware status:	FF 00000000	
Lamp counter:	17	
System test:	✓	
Filter test:	✓	
Lamp test:	✓	
Wavelength calibration:	✓	
Free internal memory space:	F3S 50176 KB	
Registered user:	Mateusz	

1.65 TIMER

This « Timer » menu allows storing in memory a waiting time before to start a measurement (Ex: Incubation time of a react ional mixture). This « Timer » acts like a stop watch.

- Press <TIMER> key
- Press F3 key [*Edit*] to enter a delay.
- Press F1 key [*Start*] to start the time count down.
- Press F2 key [*Stop*] to stop the time count down.

When the programmed delay is ended,

- The timer displays 00:00:00,
- Its status commutes from « *Running* » to « *Expired* » and
- One can hear a “bip”. Press F2 key [*Stop*] to deactivate the timer.

Timer			09/12/15 13:34
Designation	Time	Status	
User defined timer	00:00:15	Inactive	
test - 1	00:01:00	Inactive	
test - 2	00:10:00	Inactive	

Start Stop Remove Add

NOTE:

The [Add] function key is only displayed if a method is selected for which analysis timers were programmed in concentration mode. Press < timer > in the concentration mode.

1.66 SOFTWARE UPDATING

The software and method update is used to continuously update the spectrophotometer.

Remark

Only users of the user group, *Administrator* are allowed to carry out a software and method update.

The update comprises

- The newest firmware (meter software)
- New or changed method data

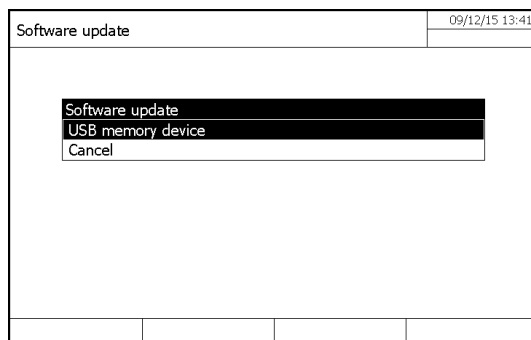
Data transfer to the spectrophotometer can be made with USB medium as a temporary storage

To realize the update, new software version or new methods have to be recording on USB medium. This one has to be connected to the spectrophotometer.

Execution

1. Connect the USB medium to the PC.
2. Unpack the contents and complete folder structure of the downloaded exe or zip file in the main directory (top level) of the USB medium (If an unpacking program such as WinZip is used, the option, "Nutz Ordnernamen" or "Use Folder Names" must be set).

3. Connect the USB medium to the spectrophotometer.
4. Switch on the spectrophotometer if necessary.

Press «▲» and «▼» keys to select « USB memory » as a source and press <START-ENTER> key.
The update process takes about five minutes. The spectrophotometer switches itself off and then on again.

Remark

If the update cannot be carried out, an error message appears on the display.

Check whether the "Update" folder with its subfolders is stored on the USB memory device (top level).

If on the photometer there is not enough free memory capacity for the update, you can create memory capacity by erasing measurement data. Save your data to a USB memory device before erasing them on the photometer.

16 MAINTENANCE

1.67 LAMP CHANGING

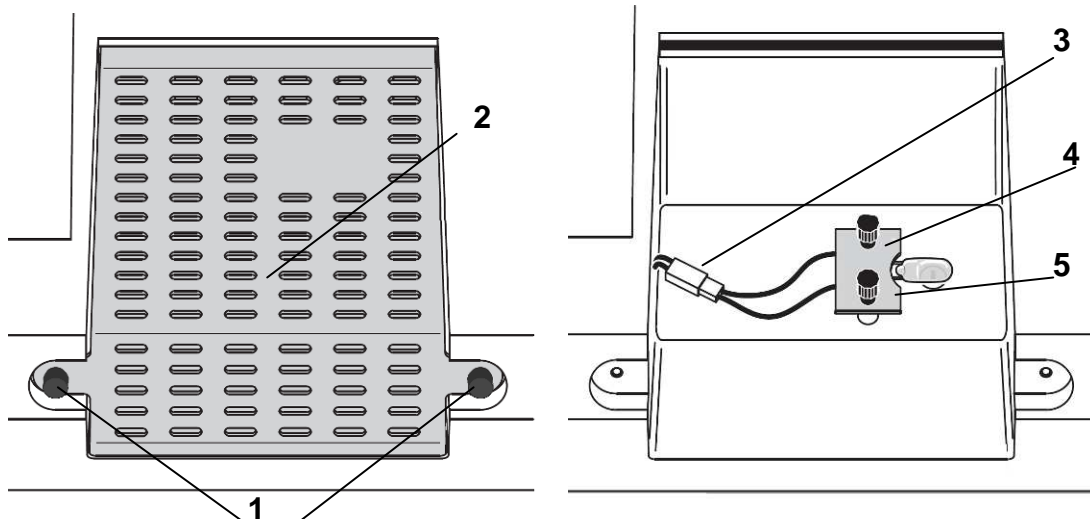
In this chapter only the change of the UviLine 9300 halogen lamp is explained. Indeed, UV-Vis UviLine spectrophotometer works with xenon lamp which time life approaches the one of the apparatus. Any time, if it prematurely stops working, then contact AQUALABO ANALYSE servicing.

The tungsten halogen lamp is a wear part with a certain average service life. It has to be replaced if defective. The spectrophotometer has a service hour counter for the lamp module.

NOTE

The replacement lamp is readily assembled as a lamp module and optically adjusted in the factory. Therefore, treat it with utmost care. Fingerprints on the glass will shorten the service life of the lamp. Do not touch the bulb of the new lamp module with bare fingers. If you have touched the bulb inadvertently, carefully clean it with a clean cloth soaked in alcohol.

The halogen lamp is behind a protection cover (Aluminum sheet) at the rear part of the spectrophotometer. To change it, proceed as follows:

WARNING

The lamp becomes very hot during operation. Do not touch the hot lamp because it can cause burns! The lamp should cool down for approx. 10 minutes before it is exchanged.

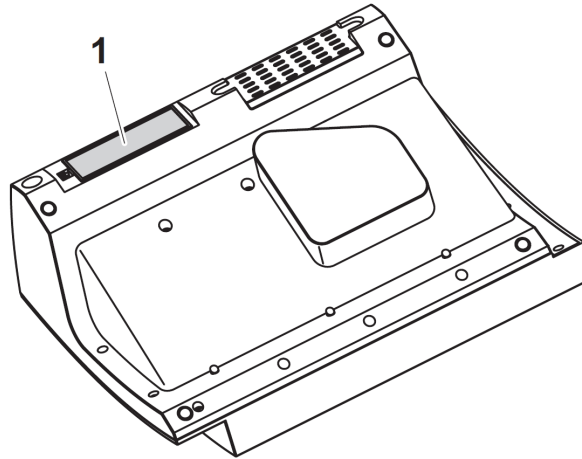
1. Switch off the spectrophotometer and disconnect it from the main power.
2. Remove the two screws (1) and remove the lamp protection cover (2).
3. Disconnect the electrical plug connection (3). Disconnecting it might take a lot of effort.
4. Unscrew the two knurled-head screws (4) and remove the defective lamp module (5).
5. Fix the new lamp module with the knurled-head screws. When doing so, the metal-plated side of the PCB must point outward, toward the knurled-head screws.
6. Reconnect the power supply cable.
7. Reinstall the protection lamp cover.
8. Reset the service hour counter for the lamp module.

1.68 BUFFER BATTERIES CHANGING

Information

Only use leak proof alkaline manganese batteries. If you leave the spectrophotometer switched on during the exchange or insert the new batteries within a minute after taking out the old ones, the date and time are retained in the spectrophotometer.

1. Turn on the spectrophotometer upside down and place it on a soft surface.
2. Remove the cover (1) of batteries compartment.
3. Remove out of the compartment the old batteries.
4. Insert the four new batteries in the battery compartment. Make sure that the poles of the batteries are in the correct position.
5. The \pm signs on the batteries must correspond to the \pm signs in the battery compartment.
6. Close the batteries compartment cover.

Batteries time life

The power consumption of the clock is very low. The lifetime of high quality batteries is at least 5 years.

1.69 CLEANING

WARNING

The housing components are made out of synthetic materials (ABS, PMMA and PC). Thus, avoid contact with acetone, ethyl alcohol and similar detergents that contain solvents. Any splashes must be wiped off immediately.

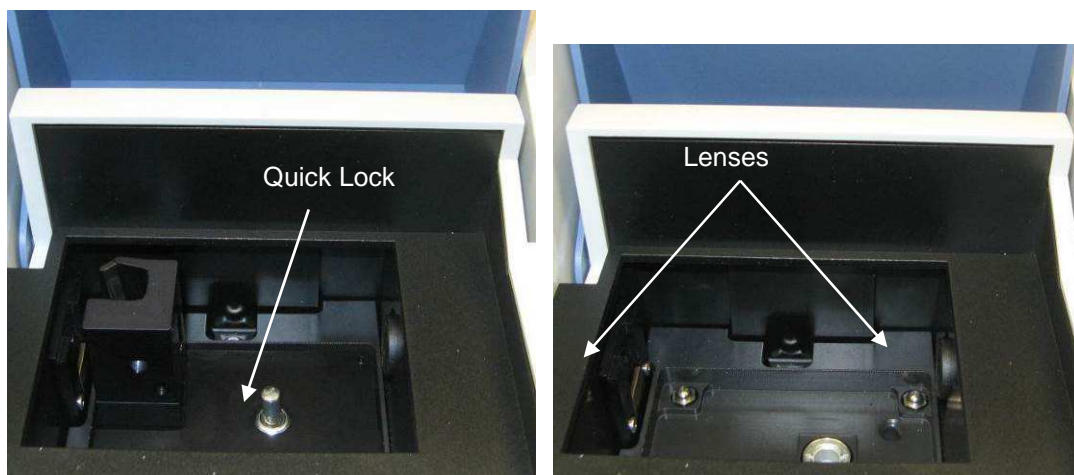
1.69.1 CLEANING YOUR SEPECTROPHOTOMETER

Clean your photometer as follows:

- If the housing surface is dirty, wipe it with a soft cloth and mild soapy water.
- Remove any chemicals splashes as soon as possible.
- For disinfection, you can use isopropanol for cleaning for a short time.

1.69.2 CLEANING THE LENSES

AQUALABO ANALYSE advises to regularly check that the 2 lenses inside the cell holder are clean. In order to access the lenses, use the Quick –Lock system to take out the accessory inserted in the cell holder (see picture below). If necessary, then proceed to the cautious cleaning of the 2 lenses using a cotton bud lightly soaked with isopropanol. Check there are no deposit/traces before use.

1.70 WHAT CAN YOU DO, IF...

1.70.1 WARNING IF CUVETTE IS BROKEN

WARNING

Cells can contain dangerous substances. If the contents are released, follow the safety instructions in the package insert. If necessary, take corresponding protective measures (protective goggles, protective gloves etc.).

CAUTION

Do not turn the photometer upside down to remove the liquid! When doing so, the liquid could come into contact with electronic components and damage the photometer.

The photometer has a drain device through which the contents of a broken cell can drain off without causing any damage.

Proceeding after a cuvette is broken:

1. Switch of the spectrophotometer and disconnect it from the main power supply.
2. Let the liquid drain off into a suitable container and dispose of it properly according to the instructions of the reagent package.
3. Carefully clean the cuvette well using a wet, lint-free cloth.
4. Let dry the cuvette well.

Remark

After recommissioning, check out the spectrophotometer for all measurements.

1.71 TROUBLESHOOTING

The spectrophotometer doesn't switch on

Cause	Remedy
- The power cable is not connected	- Connect correctly the power cable

Measurement range undercut or exceeded

Cause	Remedy
- Method not suitable.	- Select another method which suitable measuring range. - Dilute the sample.

Not stable results

Cause	Remedy
- The cover of cuvette compartment is opened	- Close the cover of the cuvette compartment
- The halogen lamp arrives at its end life	- Change the halogen lamp

Incorrect measured values

Cause	Remedy
- Cuvette dirty	- Clean the cuvette
- Wrong dilution	- Make a dilution again
- The selected method is not appropriated	- Select another method
- Erroneous zero measurement	- Perform the zero again
- Erroneous blank value	- Measure the blank again

Connected printer does not print

Cause	Remedy
- Printer not suitable	- Connect a printer that can interpret the printer control language PCL-3

Data transmission to USB memory device does not work

Cause	Remedy
- Connected USB memory device was not recognized - The USB memory device has been formatted to a file system which is not supported, e. g. (NTFS)	- Use other USB memory device - Format the USB memory device to the FAT 32 file system

Incorrect measured values

Cause	Remedy
- Cuvette dirty	- Clean the cuvette
- Wrong dilution	- Make a dilution again
- The selected method is not appropriated	- Select another method
- Erroneous zero measurement	- Perform the zero again
- Erroneous blank value	- Measure the blank again

17 ACCESSORIES

List of manual and automatic accessories available on UviLine instruments

1.72 10MM CELL HOLDER (REF. 70VI0604)

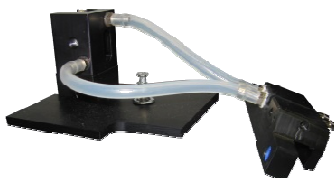
Single cell holder up to 10mm optical path length. The 10mm cell holder guarantees a perfect positioning for all types of cells. It can be used with micro-cells up to 50 μ l. Quick and easy installation thanks to the Quick-Lock system.

1.73 PORTE CUVE JUSQU'A 100MM (REF. 70VI0605)

Tank holder up to 100mm of optical thickness for tanks (comes as standard with PASTEL UviLine devices). Suitable for 5/10/20/50/100mm tanks and 16mm tubes, it rigorously covers all manual uses of the UviLines that require high reading sensitivity. Quick and easy installation due to the Quick-Lock system.

1.74 PORTE CUVE 10MM THERMOSTATABLE (REF. 70VI0602)


Thermostatable 10mm cell holder with integrated thermostatable function. Allows all standard cell with a 10mm optical path length. Quick and easy installation thanks to the Quick-Lock system. Thermostatable by external bath.

1.75 PORTE TUBE 16MM (REF. 70VI0607)

16mm tube holder. Can be used in combination with micro-methods and especially designed for the use of 16mm diameters tubes. Quick and easy installation thanks to the Quick-Lock system and the sketch indicating how to position the tube.

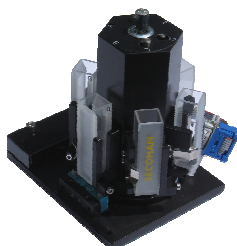
1.76 AUTOMATIC CELL CHANGER TURRET (REF. 70VI0600)

Designed to analyze 5 samples and 1 blank, the 5+1 turret find its place in automation applications and guarantees measurements stability with unique advantages:

- Removable and easy to re-insert turret for convenient cells set up
- Accurate positioning for low volume cells

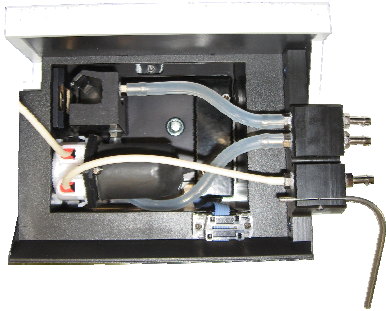
1.77 SIPPER (REF. 70VI0601)

Compact, the sipper is delivered standard with a peristaltic pump integrated to the cell holder allowing:

- Programmable sucking from 500 μ l to 3000 μ l
- Use of flowthrough cells from 30 μ l to 450 μ l
- It makes manipulation safer and higher in productivity

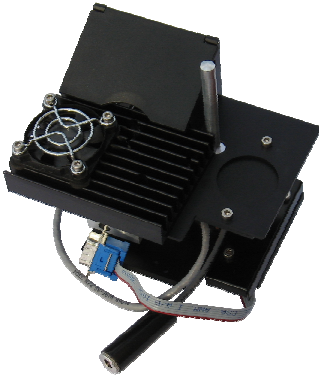1.78 EFFET PELTIER (REF. 70VI0603)

- T°C set up between 10°C and 60°C
- Accuracy: \pm 0.5°C
- Sipper option
- Fast, ultra compact and fully instrument driven

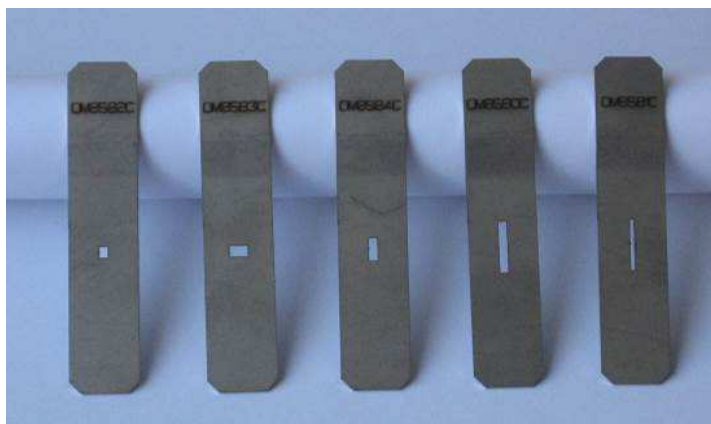
1.79 SIPPER WITH PELTIER EFFECT (REF. 70VI0610)

Compact, this accessory is equipped with a peristaltic pump cassette integrated to the cell holder and a Peltier module temperature regulating system. Temperature programming between 10°C and 60°C

- Accuracy: + 0.5°C
- Aspiration programming: from 250 μ l to 2000 μ l
- Cells: from 30 μ l to 450 μ l
- It secures and increases productivity handling
- Very compact, fast and completely controlled by the device

1.80 DIAPHRAGMS FOR MICRO-CELLS (REF. 70VI0608)

Reference: 70VI0608 (optional accessory)
Designation: Batch of 5 diaphragms for micro-cell

Reference 70VI0608 includes the following 5 diaphragms: OM8582 : size 1,8 x 2,3 mm ; OM8583 : size 3,8 x 2,3 mm ; OM8584 : size 1,8 x 4,8 mm ; OM8580 : size 1,8 x 11 mm ; OM8581 : size 0,9 x 11 mm

Use: these diaphragms are to be used with micro-cells without black walls. They help concentrate the beam on the sample through the hole in the metal strip, without the light being refracted in all directions.

You will find here below the list of SECOMAM cells for which we advise to use the diaphragms. This list indicates the reference of the diaphragm to be used with the micro-cells.

Cell reference	Diaphragm reference
0GQ53Z0	OM8580
0GG53Z0	OM8580
0GQ54Z0	OM8580
0GG54Z0	OM8580
0GQ123Z0	OM8581
0GG123Z0	OM8581
0GQ124Z0	OM8581
0GG124Z0	OM8581
0GQ133Z0	OM8580
0GG133Z0	OM8580
0GQ134Z0	OM8580
0GG134Z0	OM8580
0GQ233Z0	OM8581
0GG233Z0	OM8581
0GQ234Z0	OM8581
0GG234Z0	OM8581
0GQ243Z0	OM8580
0GG243Z0	OM8580

Cell reference	Diaphragm reference
0GQ244Z0	OM8580
0GG244Z0	OM8580
0GQ253Z0	OM8580
0GG253Z0	OM8580
0GQ254Z0	OM8580
0GG254Z0	OM8580
0GQ283Z0	OM8580
0GG283Z0	OM8580
0GQ284Z0	OM8580
0GG284Z0	OM8580
0GQ373Z0	OM8581
0GG373Z0	OM8581
0GQ374Z0	OM8581
0GG374Z0	OM8581
0GQ383Z0	OM8580
0GG383Z0	OM8580
0GQ384Z0	OM8580
0GG384Z0	OM8580

Cell reference	Diaphragm reference
0GQ424Z0	OM8581
0GG424Z0	OM8581
0GQ434Z0	OM8580
0GG434Z0	OM8580
0GQ444Z0	OM8580
0GG444Z0	OM8580
0GQ464Z0	OM8581
0GG464Z0	OM8581
0GQ474Z0	OM8580
0GG474Z0	OM8580
0GQ484Z0	OM8580
0GG484Z0	OM8580
0GQ634Z8.5	OM8582
0GG634Z15	OM8582
0GQ634Z8.5	OM8582
0GG634Z15	OM8582

Cell reference	Diaphragm reference
0GQ644Z8.5	OM8583
0GG644Z15	OM8583
0GQ644Z8.5	OM8583
0GG644Z15	OM8583
0GQ653Z8.5	OM8582
0GG653Z15	OM8582
0GQ653Z8.5	OM8582
0GG653Z15	OM8582
0GQ654Z8.5	OM8583
0GG654Z15	OM8583
0GQ654Z8.5	OM8583
0GG654Z15	OM8583
0GQ674Z8.5	OM8582
0GG674Z15	OM8582
0GQ674Z15	OM8582
0GG674Z8.5	OM8582
0GG674Z15	OM8582

SECOMAM also sells a large range of cells for which you will find here below the list of our best-sellers.
For our full list of available cells, contact us on analytics.info-fr@xyleminc.com

Our Best-sellers :

References	Description (packaging: one cell)
0GQ4Z0	Standard quartz cell with lid 10mm 3,5ml 45*12,5*12,5
0GG4Z0	Standard glass cell with lid 10mm 3,5ml 45*12,5*12,5
0GQ5Z0	Standard quartz cell with lid 20mm 7ml 45*12,5*22,5
0GG5Z0	Standard glass cell with lid 20mm 7ml 45*12,5*22,5
0GQ8Z0	Standard quartz cell with lid 50 mm 17,5ml 45*12,5*52,5
0GG8Z0	Standard glass cell with lid 50 mm 17,5ml 45*12,5*52,5
0GQ14Z0	Standard quartz cell with teflon lid 10mm 3,5ml 45*12,5*12,5
0GG14Z0	Standard glass cell with teflon lid 10mm 3,5ml 45*12,5*12,5
0GQ15Z0	Standard quartz cell with teflon lid 20mm 7ml 45*12,5*22,5
0GG15Z0	Standard glass cell with teflon lid 20mm 7ml 45*12,5*22,5
0GQ18Z0	Standard quartz cell with teflon lid 50 mm 17,5ml 45*12,5*52,5
0GG18Z0	Standard glass cell with teflon lid 50 mm 17,5ml 45*12,5*52,5
0GQ74Z8.5	Glass flowthrough cell 10 mm 0,48ml 45*12,5*12,5
0GQ74Z15	Glass flowthrough cell 10 mm 0,48ml 45*12,5*12,5
0GG74Z8.5	Glass flowthrough cell 10 mm 0,48ml 45*12,5*12,5
0GG74Z15	Glass flowthrough cell 10 mm 0,48ml 45*12,5*12,5
0GQ244Z0	Micro quartz cell with frosted walls and with lid 10 mm 0,7ml 45*12,5*12,5
0GG244Z0	Micro glass cell with frosted walls and with lid 10 mm 0,7ml 45*12,5*12,5
0GQ245Z0	Micro quartz cell with frosted walls and with lid 20 mm 1,4ml 45*12,5*22,5
0GG245Z0	Micro glass cell with frosted walls and with lid 20 mm 1,4ml 45*12,5*22,5
0GQ254Z0	Micro quartz cell with frosted walls and with lid 10 mm 1ml 45*12,5*12,5
0GG254Z0	Micro glass cell with frosted walls and with lid 10 mm 1ml 45*12,5*12,5
0GQ255Z0	Micro quartz cell with frosted walls and with lid 20 mm 2,1ml 45*12,5*22,5
0GG255Z0	Micro glass cell with frosted walls and with lid 20 mm 2,1ml 45*12,5*22,5
0GQ264Z0	Semi micro quartz cell with frosted walls and with lid 10 mm 1,4ml 45*12,5*12,5
0GG264Z0	Semi micro glass cell with frosted walls and with lid 10 mm 1,4ml 45*12,5*12,5
0GQ265Z0	Semi micro quartz cell with frosted walls and with lid 20 mm 2,8ml 45*12,5*22,5
0GG265Z0	Semi micro glass cell with frosted walls and with lid 20 mm 2,8ml 45*12,5*22,5
0GQ268Z0	Semi micro quartz cell with frosted walls and with lid 50 mm 7ml 45*12,5*52,5
0GG268Z0	Semi micro glass cell with frosted walls and with lid 50 mm 7ml 45*12,5*52,5
0GQ653Z8.5	Quartz micro cell with black walls 10 mm 0,05 ml 13*12,5*12,5
0GQ653Z15	Quartz micro cell with black walls 10 mm 0,05 ml 13*12,5*12,5
0GG653Z8.5	Glass micro cell with black walls 10 mm 0,05 ml 13*12,5*12,5
0GG653Z15	Glass micro cell with black walls 10 mm 0,05 ml 13*12,5*12,5

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The diagram illustrates the navigation flow between different screens of the UviLine device. It starts with the 'Home' screen, which shows five main modes: Concentration, Absorbance /% transmission, Multi-longueurs d'onde, Spectre, and Cinétique. A central navigation pad with four arrow keys and ESC/START ENTER buttons is shown. Arrows indicate the following paths:

- From the 'Home' screen, pressing F2 leads to the 'Configuration' screen.
- From the 'Home' screen, pressing F1 leads to the 'Concentration' screen.
- From the 'Home' screen, pressing F3 leads to the 'Absorbance' screen.
- From the 'Home' screen, pressing F4 leads to the 'Spectre' screen.

Home Screen (01/01/15 12:01): Shows icons for Concentration (C), Absorbance (% transmission) (λ), Multi-longueurs d'onde (Σ), Spectre (λ), and Cinétique (t). Buttons for Configuration, Info, F1, F2, F3, and F4 are visible.

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Concentration Screen (04/16/15 14:40): Displays a large reading of 12.24 mg/l and Absorbance 0.562. It also shows '1000: Nitrate' and 'NO3 10.00 - 30.00 mg/l'. Buttons for Paramètre and Liste de méth. are at the bottom.

Absorbance Screen (04/16/15 15:12): Shows a reading of 0.079 and 83.3%. A table of data is displayed:

0.079	83.5 %
0.081	82.9 %
0.080	83.2 %
0.078	83.6 %

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Spectre Screen (04/16/15 15:52): Displays a graph of Absorbance vs. Longueur d'onde [nm]. The x-axis ranges from 400 to 600 nm. The y-axis ranges from 0.0 to 0.6. A peak is visible at 537.0 nm with an absorbance of 0.886. Buttons for Paramètre, Calculs, Zoom, and Ouvrir are at the bottom.

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
Edit method		04/16/15 14:41
Number	1002	
Wavelength	190 nm	
Resoluti	Calibration curve	0.01
Unit	Measure standard solutions	mg/l
Citation	Enter formula	
Designa	Enter value pairs	
Version		
Calibration curve	Measure standard solutions	
Cell changer mode	Single cell	
Back		Next

Edit method		04/16/15 14:43
Blank value		0.010
Std. 1	10.00 mg/l	0.198
Std. 2	20.00 mg/l	0.405
Std. 3	30.00 mg/l	0.810
Back		Next

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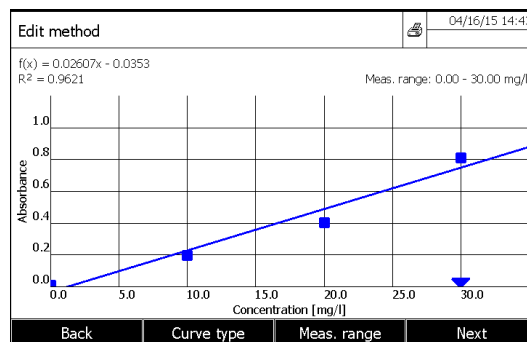
Edit method		04/16/15 14:44
Number	1002	
Wavelength	190 nm	
Resoluti	Calibration curve	0.01
Unit	Measure standard solutions	mg/l
Citation	Enter formula	
Designa	Enter value pairs	
Version		
Calibration curve	Enter value pairs	
Cell changer mode	Single cell	
Back		Next


Edit method		04/16/15 14:46
$c = a_0 + a_1 \cdot A + a_2 \cdot A^2 + a_3 \cdot A^3 + a_4 \cdot A^4 + a_5 \cdot A^5$		
a0		
a1		
a2		
a3		
a4		
a5		
Lower limit of measuring range		
Upper limit of measuring range		
Back		Next

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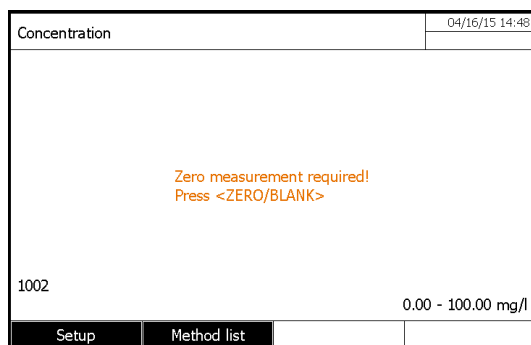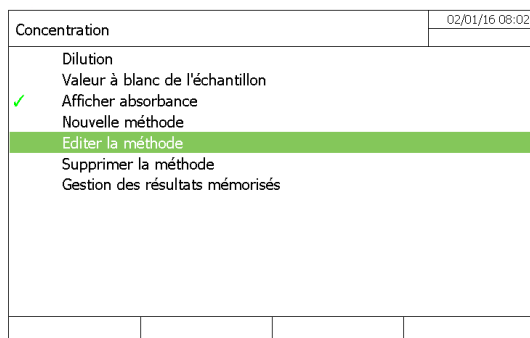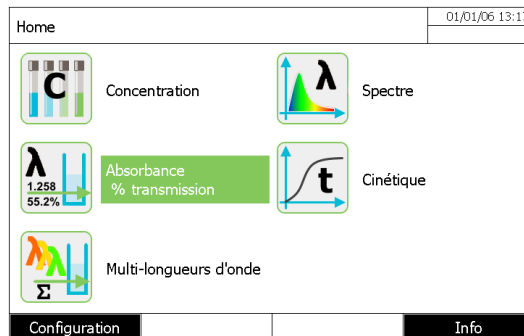
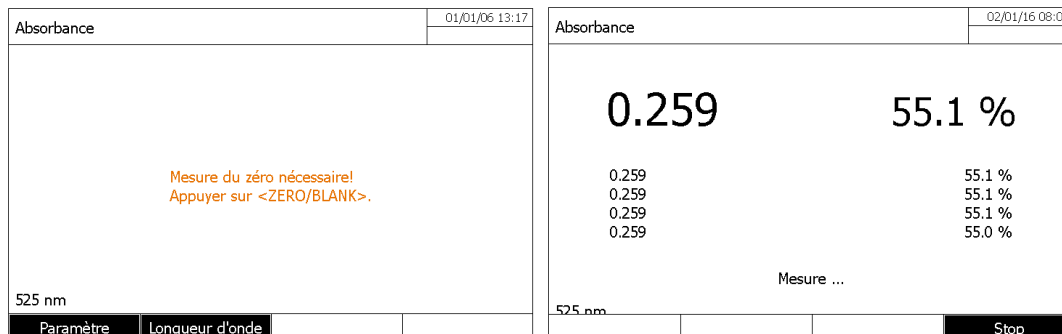
```

UviLine 9300 7449001 1.14
Administrator

Time of measurement:      07/01/2015 14:05:46
Wavelength:              525 nm
Measured value:          0.278  52.8 %
    
```

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© 2015 AQUALABO-GROUP - All rights reserved